Operating Instructions Memosens CLS82E

Hygienic conductivity sensor Digital with Memosens technology







Table of contents Memosens CLS82E

Table of contents

1 1.1 1.2 1.3	About this document3Warnings3Symbols3Documentation3
2 2.1 2.2 2.3 2.4 2.5	Basic safety instructions4Requirements for the personnel4Intended use4Workplace safety4Operational safety5Product safety5
3	Incoming acceptance and
3.1 3.2 3.3	product identification5Incoming acceptance5Product identification6Scope of delivery6
4 4.1 4.2	Mounting7Mounting requirements7Post-mounting check10
5 5.1 5.2 5.3	Electrical connection10Connecting the sensor11Ensuring the degree of protection11Post-connection check11
6	Commissioning 12
7 7.1 7.2	Maintenance12Cleaning the sensor12Sensor calibration13
8 8.1 8.2 8.3 8.4	Repair 13 General notes 13 Spare parts 14 Return 14 Disposal 14
9 9.1 9.2	Accessories15Measuring cable15Calibration solutions15
10 10.1 10.2	Technical data15Input15Performance characteristics16

Index 19			
10.5	Mechanical construction	18	
10.4	Process	17	
10.3	Environment	17	

Memosens CLS82E About this document

1 About this document

1.1 Warnings

Structure of information	Meaning
▲ DANGER Causes (/consequences) If necessary, Consequences of non- compliance (if applicable) Corrective action	This symbol alerts you to a dangerous situation. Failure to avoid the dangerous situation will result in a fatal or serious injury.
▲ WARNING Causes (/consequences) If necessary, Consequences of non- compliance (if applicable) Corrective action	This symbol alerts you to a dangerous situation. Failure to avoid the dangerous situation can result in a fatal or serious injury.
Causes (/consequences) If necessary, Consequences of non- compliance (if applicable) Corrective action	This symbol alerts you to a dangerous situation. Failure to avoid this situation can result in minor or more serious injuries.
NOTICE Cause/situation If necessary, Consequences of non- compliance (if applicable) Action/note	This symbol alerts you to situations which may result in damage to property.

1.2 Symbols

Additional information, tips
Permitted or recommended
Not permitted or not recommended
Reference to device documentation
Reference to page
Reference to graphic
Result of a step

1.3 Documentation



 $Technical\ Information\ Memosens\ CLS82E,\ TIO1529C$



Special Documentation for hygienic applications, SD02751C

In addition to these Operating Instructions, an XA with "Safety instructions for electrical apparatus in the hazardous area" is also included with sensors for use in the hazardous area.


▶ Please follow instructions on use in the hazardous area carefully.

Basic safety instructions Memosens CLS82E

2 Basic safety instructions

2.1 Requirements for the personnel

- Installation, commissioning, operation and maintenance of the measuring system may be carried out only by specially trained technical personnel.
- The technical personnel must be authorized by the plant operator to carry out the specified activities.
- The electrical connection may be performed only by an electrical technician.
- The technical personnel must have read and understood these Operating Instructions and must follow the instructions contained therein.
- Faults at the measuring point may only be rectified by authorized and specially trained personnel.

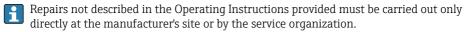

2.2 Intended use

The Memosens CLS82E conductivity sensor is used to measure low to high conductivity of liquids in applications with hygienic requirements.

The broad measuring range means the device can be used in a large number of applications, e.g.:

- Phase separation of water/product mixtures
- Phase separation of product/product mixtures
- Monitoring of rinsing processes
- Fermentations
- Monitoring of water bodies
- Concentration measurement of bases and acids (consider the material resistance properties!)
- Monitoring product quality

The digital sensor is used with the Liquiline CM44x or Liquiline CM42.

Use of the device for any purpose other than that described, poses a threat to the safety of people and of the entire measuring system and is therefore not permitted.

The manufacturer is not liable for damage caused by improper or non-designated use.

2.3 Workplace safety

As the user, you are responsible for complying with the following safety conditions:

- Installation guidelines
- Local standards and regulations
- ullet Regulations for explosion protection

Electromagnetic compatibility

- The product has been tested for electromagnetic compatibility in accordance with the applicable international standards for industrial applications.
- The electromagnetic compatibility indicated applies only to a product that has been connected in accordance with these Operating Instructions.

2.4 Operational safety

Before commissioning the entire measuring point:

- 1. Verify that all connections are correct.
- 2. Ensure that electrical cables and hose connections are undamaged.
- 3. Do not operate damaged products, and protect them against unintentional operation.
- 4. Label damaged products as defective.

During operation:

► If faults cannot be rectified: products must be taken out of service and protected against unintentional operation.

2.5 Product safety

The product is designed to meet state-of-the-art safety requirements, has been tested, and left the factory in a condition in which it is safe to operate. The relevant regulations and international standards have been observed.

3 Incoming acceptance and product identification

3.1 Incoming acceptance

- 1. Verify that the packaging is undamaged.
 - Wotify the supplier of any damage to the packaging.
 Keep the damaged packaging until the issue has been resolved.
- 2. Verify that the contents are undamaged.
 - Notify the supplier of any damage to the delivery contents. Keep the damaged goods until the issue has been resolved.
- 3. Check that the delivery is complete and nothing is missing.
 - ightharpoonup Compare the shipping documents with your order.
- **4.** Pack the product for storage and transportation in such a way that it is protected against impact and moisture.
 - The original packaging offers the best protection.

 Make sure to comply with the permitted ambient conditions.

If you have any questions, please contact your supplier or your local Sales Center.

3.2 Product identification

3.2.1 Nameplate

The nameplate provides you with the following information on your device:

- Manufacturer identification
- Extended order code
- Serial number
- Safety information and warnings
- ► Compare the information on the nameplate with the order.

3.2.2 Product identification

Product page

www.endress.com/cls82e

Interpreting the order code

The order code and serial number of your product can be found in the following locations:

- On the nameplate
- In the delivery papers

Obtaining information on the product

- 1. Open www.endress.com.
- 2. Call up the site search (magnifying glass).
- 3. Enter a valid serial number.
- 4. Search.
 - ► The product structure is displayed in a popup window.
- 5. Click on the product image in the popup window.
 - A new window (**Device Viewer**) opens. All of the information relating to your device is displayed in this window as well as the product documentation.

Manufacturer address

Endress+Hauser Conducta GmbH+Co. KG Dieselstraße 24 D-70839 Gerlingen

3.3 Scope of delivery

The scope of delivery includes:

- Sensor in the version ordered
- Operating Instructions

Memosens CLS82E Mounting

4 Mounting

4.1 Mounting requirements

4.1.1 Hygiene-compliant installation


► The use of an EHEDG-certified assembly is a prerequisite for the easy-to-clean installation of a 12-mm sensor in accordance with EHEDG requirements.

- ► Furthermore, the instructions regarding the hygienic installation and operation of the assembly in the relevant Operating Instructions must be adhered to.
- Easily cleanable installation of equipment according to the criteria of the EHEDG must be free of dead legs.
- ▶ If a dead leg is unavoidable, it shall be kept as short as possible. Under no circumstances shall the length of a dead leg L exceed the pipe's inner diameter D minus the equipment's enveloping diameter d. The condition $L \le D d$ applies.
- ► Furthermore, the dead leg must be self-draining, so neither product nor process fluids are retained therein.
- Within tank installations, the cleaning device must be located so that it directly flushes the dead leq.
- ► For further reference, see the recommendations concerning hygienic seals and installations in EHEDG Doc. 10 and the Position Paper: "Easy cleanable Pipe couplings and Process connections".

For 3-A-compliant installation, please observe the following:

- ▶ After the device has been mounted, hygienic integrity must be guaranteed.
- ► 3-A-compliant process connections must be used.

4.1.2 Installation factors for assemblies

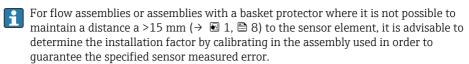


▶ Prior to installation:

Remove the black protective cap from the sensor element.

Symmetrical installation is recommended in order to guarantee linearity. The distance to the side walls and opposite walls must be at least 15 mm.

Mounting Memosens CLS82E

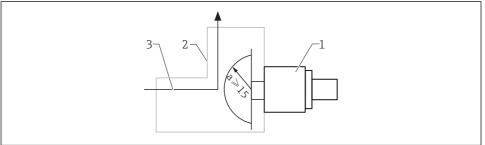

40024421

 \blacksquare 1 Minimum distance between pipe and end of the measuring cell


- 1 Sensor
- 2 Pipe
- 3 Direction of flow

The ionic current in the liquid is affected by the walls in confined installation conditions. This effect is compensated by what is referred to as the installation factor. The installation factor can be entered in the transmitter for the measurement or the cell constant is corrected by multiplying by the installation factor.

The value of the installation factor depends on the diameter and the conductivity of the pipe nozzle as well as the sensor's distance to the wall. The installation factor can be disregarded (f = 1.00) if the distance to the wall is sufficient (a > 15 mm). If the distance to the wall is smaller, the installation factor increases for electrically insulating pipes (f > 1) and decreases for electrically conductive pipes (f < 1). The installation factor can be determined using calibration solutions

- ► Ensure that the electrodes are fully immersed in the medium during measurement. Ideally, medium should flow to the measuring cell from the front.
 - Any other installation position can cause air pockets to occur or the buildup of solid impurities.


Memosens CLS82E Mounting

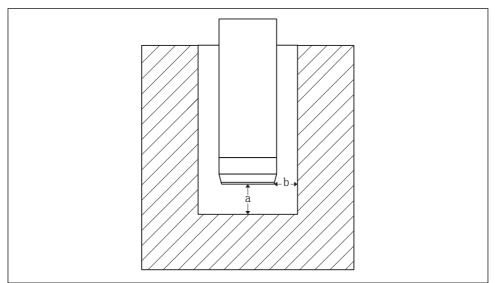

A0024626

■ 2 Schematic drawing of the sensor in confined installation conditions


- a Wall distance
- b Gap width

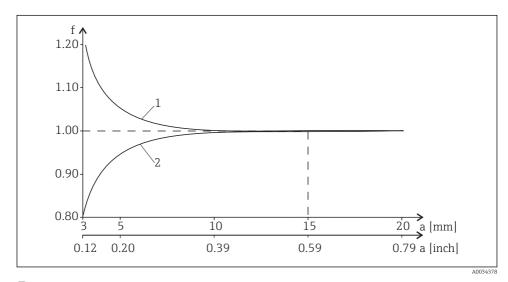

 \blacksquare 3 Relationship between installation factor f and wall distance a

- 1 Electrically insulating pipe wall
- 2 Electrically conductive pipe wall

Electrical connection Memosens CLS82E

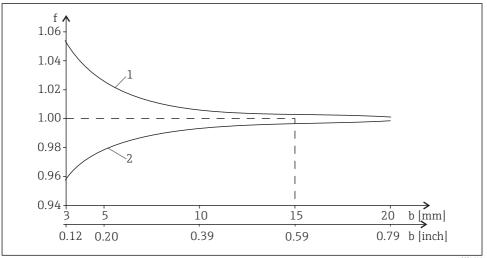

A0024616

 \blacksquare 4 Relationship between installation factor f and gap width b

- 1 Electrically insulating pipe wall
- 2 Electrically conductive pipe wall

4.2 Post-mounting check

- 1. Are the sensor and cable undamaged?
- 2. Is the sensor installed in the process connection and is not suspended from the cable?

5 Electrical connection

A WARNING

Device is live!


Incorrect connection may result in injury or death!

- ► The electrical connection may be performed only by an electrical technician.
- ► The electrical technician must have read and understood these Operating Instructions and must follow the instructions contained therein.
- ▶ **Prior** to commencing connection work, ensure that no voltage is present on any cable.

Memosens CLS82E Electrical connection

5.1 Connecting the sensor

The electrical connection of the sensor to the transmitter is established using the measuring cable CYK10.

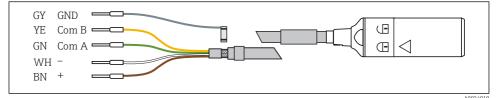

M : 11 CW//10

■ 5 Measuring cable CYK10

5.2 Ensuring the degree of protection

Only the mechanical and electrical connections which are described in these instructions and which are necessary for the required, designated use, may be carried out on the device delivered.

► Exercise care when carrying out the work.

Otherwise, the individual types of protection (Ingress Protection (IP), electrical safety, EMC interference immunity) agreed for this product can no longer be guaranteed due, for example to covers being left off or cable (ends) that are loose or insufficiently secured.

5.3 Post-connection check

WARNING

Connection errors

The safety of people and of the measuring point is at risk! The manufacturer does not accept any responsibility for errors that result from failure to comply with the instructions in this manual.

Put the measuring point into operation only if you can answer yes to all the following questions.

Product status and specifications

▶ Are the sensor and cable free from damage on the outside?

Electrical connection


- ▶ Is the installed cable strain-relieved and not twisted?
- ► Is a sufficient length of the cable cores stripped, and are the cores correctly positioned in the terminal on the transmitter?
- ► Are all plug-in terminals on the transmitter securely engaged?
- ► Are all cable entries mounted on the transmitter, tightened and leak-tight?

Commissioning Memosens CLS82E

6 Commissioning

Prior to initial commissioning, ensure that:

- The sensor is correctly installed
- The electrical connection is correct
- 1. Check the temperature compensation and damping settings on the transmitter.

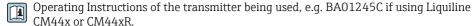

WARNING

Escaping process medium

Risk of injury from high pressure, high temperatures or chemical hazards!

- ► Before applying pressure to an assembly with cleaning system, ensure that the system has been connected correctly.
- ► If you cannot reliably establish the correct connection, do not install the assembly in the process.

If using an assembly with automatic cleaning function:

- 2. Check that the cleaning medium (water or air, for example) is connected correctly.
- 3. Following commissioning:

Maintain the sensor at regular intervals.

- This is the only way to ensure reliable measurements.
- As the sensor can be operated with a nominal pressure greater than 1 bar (15 psi), it has been registered according to CSA B51 ("Boiler, pressure vessel, and pressure piping code"; category F) with a CRN (Canadian Registration Number) in all Canadian provinces.

The CRN is located on the nameplate.

7 Maintenance

7.1 Cleaning the sensor

A CAUTION

Corrosive chemicals

Risk of chemical burns to the eyes and skin and risk of damage to clothing and equipment!

- ► It is absolutely essential to protect the eyes and hands properly when working with acids, alkalis and organic solvents!
- ► Wear protective goggles and safety gloves.
- $\,\blacktriangleright\,$ Clean away splashes on clothes and other objects to prevent any damage.
- ► Comply with instructions in the safety data sheets for the chemicals used.

Memosens CLS82E Repair

A WARNING

Thiocarbamide

Harmful if swallowed! Limited evidence of carcinogenicity! Possible risk of harm to the unborn child! Dangerous for the environment with long-term effects!

- ▶ Wear protective goggles, protective gloves and appropriate protective clothing.
- ▶ Avoid all contact with the eyes, mouth and skin.
- ▶ Avoid discharge into the environment.

Clean away fouling on the sensor as follows depending on the type of fouling:

- 1. Oily and greasy films:
 - Clean with fat solvent, e.g. alcohol, or hot water and agents containing surfactants (alkaline) (e.g. dishwashing detergent).
- 2. Lime and metal hydroxide buildup and low solubility (lyophobic) organic buildup: Dissolve buildup with diluted hydrochloric acid (3 %) and then rinse thoroughly with plenty of clear water.
- 3. Sulfidic buildup (from flue gas desulfurization or wastewater treatment plants):
 Use a mixture of hydrochloric acid (3 %) and thiocarbamide (commercially available) and then rinse thoroughly with plenty of clear water.
- 4. Buildup containing proteins (e.g. food industry):

 Use a mixture of hydrochloric acid (0.5 %) and pepsin (commercially available) and then rinse thoroughly with plenty of clear water.
- 5. Readily soluble biological buildup: Rinse with pressurized water.

After cleaning, rinse the sensor thoroughly with water.

7.2 Sensor calibration

Wall distance:

When calibrating, ensure that there is a minimum distance of 15 mm to the base and walls of the calibration vessel.

8 Repair

8.1 General notes

The repair and conversion concept provides for the following:

- The product has a modular design
- Spare parts are grouped into kits which include the associated kit instructions
- Only use original spare parts from the manufacturer

Repair Memosens CLS82E

- Repairs are carried out by the manufacturer's Service Department or by trained users
- Certified devices can only be converted to other certified device versions by the manufacturer's Service Department or at the factory
- Observe applicable standards, national regulations, Ex documentation (XA) and certificates
- 1. Carry out the repair according to the kit instructions.
- 2. Document the repair and conversion and enter, or have entered, in the Life Cycle Management tool (W@M).

8.2 Spare parts

Device spare parts that are currently available for delivery can be found on the website:

www.endress.com/device-viewer

▶ Quote the serial number of the device when ordering spare parts.

8.3 Return

The product must be returned if repairs or a factory calibration are required, or if the wrong product was ordered or delivered. As an ISO-certified company and also due to legal regulations, Endress+Hauser is obliged to follow certain procedures when handling any returned products that have been in contact with medium.

To ensure the swift, safe and professional return of the device:

▶ Refer to the website www.endress.com/support/return-material for information on the procedure and conditions for returning devices.

8.4 Disposal



If required by the Directive 2012/19/EU on waste electrical and electronic equipment (WEEE), the product is marked with the depicted symbol in order to minimize the disposal of WEEE as unsorted municipal waste. Do not dispose of products bearing this marking as unsorted municipal waste. Instead, return them to Endress+Hauser for disposal under the applicable conditions.

Memosens CLS82E Accessories

9 Accessories

The following are the most important accessories available at the time this documentation was issued.

► For accessories not listed here, please contact your Service or Sales Center.

9.1 Measuring cable

Memosens data cable CYK10

- For digital sensors with Memosens technology
- Product Configurator on the product page: www.endress.com/cyk10



Technical Information TI00118C

Memosens data cable CYK11

- Extension cable for digital sensors with Memosens protocol
- Product Configurator on the product page: www.endress.com/cyk11



Technical Information TI00118C

9.2 Calibration solutions

Conductivity calibration solutions CLY11

Precision solutions referenced to SRM (Standard Reference Material) by NIST for qualified calibration of conductivity measuring systems in accordance with ISO 9000

- CLY11-A, 74 μS/cm (reference temperature 25 °C (77 °F)), 500 ml (16.9 fl.oz)
 Order No. 50081902
- CLY11-B, 149.6 μ S/cm (reference temperature 25 °C (77 °F)), 500 ml (16.9 fl.oz) Order No. 50081903
- CLY11-C, 1.406 mS/cm (reference temperature 25 °C (77 °F)), 500 ml (16.9 fl.oz)
 Order No. 50081904
- CLY11-D, 12.64 mS/cm (reference temperature 25 °C (77 °F)), 500 ml (16.9 fl.oz)
 Order No. 50081905
- CLY11-E, 107.00 mS/cm (reference temperature 25 °C (77 °F)), 500 ml (16.9 fl.oz)
 Order No. 50081906



Technical Information TI00162C

10 Technical data

10.1 Input

10.1.1 Measured variables

- Conductivity
- Temperature

Technical data Memosens CLS82E

10.1.2 Measuring ranges

Conductivity 1)

 $1 \mu S/cm$ to 500 mS/cm

1) In relation to water at 25 °C (77 °F)

Temperature

-5 to 140 °C (23 to 284 °F)

10.1.3 Cell constant

 $k = 0.57 \text{ cm}^{-1}$

10.1.4 Temperature compensation

Pt1000 (Class A according to IEC 60751)

10.2 Performance characteristics

10.2.1 Measuring uncertainty

Each individual sensor is factory-measured in a solution with approx. $50 \,\mu\text{S/cm}$ using a reference measuring system traceable to NIST or PTB. The exact cell constant is entered into the manufacturer inspection certificate supplied. The uncertainty of measurement in determining the cell constant is $1.0 \,\%$.

10.2.2 Response time

Conductivity

 $t_{95} \le 2 s$

Temperature 1)

With Pg 13.5 or Clamp $t_{90} \le 16 \text{ s}^{2)}$ With other process connection $t_{90} \le 28 \text{ s}^{2)}$

- 1) DIN VDI/VDE 3522-2 (0.3 m/s laminar)
- 2) With temperature prediction activated as standard

10.2.3 Maximum measured error

Conductivity

In the range 1 μ S/cm to 1 mS/cm $^{1)}$ \leq 2 % of reading In the range 1 mS/cm to 500 mS/cm $^{1)}$ \leq 4 % of reading

Temperature

With Pg 13.5 or Clamp \leq 0.5 K, in measuring range -5 to 100 $^{\circ}$ C (23 to

212 °F)

 \leq 1.0 K, in measuring range 100 to 140 °C (212 to

284°F)

With other process connection ≤ 1.0 K, in measuring range -5 to 140 °C (23 to

284°F)

1) In as-delivered state (factory adjustment at 50 μ S/cm)

Memosens CLS82E Technical data

10.2.4 Repeatability

Conductivity $\leq 0.2 \%$ of reading, in specified measuring range

Temperature $\leq 0.05 \text{ K}$

10.3 Environment

10.3.1 Ambient temperature

-20 to 60 °C (-4 to 140 °F)

10.3.2 Storage temperature

-25 to +80 °C (-10 to +180 °F)

10.3.3 Humidity

5 to 95 %

10.3.4 Degree of protection

IP 68 / NEMA type 6P (1.9 m water column, 20 °C, 24 h)

10.4 Process

10.4.1 Process temperature

Normal operation: $-5 \text{ to } 120 \,^{\circ}\text{C} (23 \text{ to } 248 \,^{\circ}\text{F})$


Sterilization (max. 45 min.): Max. $140 \,^{\circ}\text{C}$ (284 $^{\circ}\text{F}$) at 6 bar (87 psi)

10.4.2 Process pressure

17 bar (247 psi) at 20 °C (68 °F) 9 bar (131 psi) at 120 °C (248 °F)

Technical data Memosens CLS82E

10.4.3 Temperature/pressure ratings

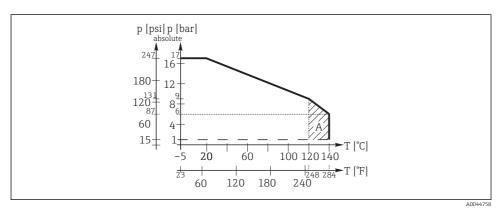

■ 6 Pressure/temperature ratings

A Can be sterilized for a short time (45 min.)

10.5 Mechanical construction

10.5.1 Weight

Depending on the version, e.g.

Process connection Pg 13.5: 0.06 to 0.09 kg (0.13 to 0.20 lbs)

• Process connection G1 or NPT: approx. 0.9 kg (1.98 lbs)

10.5.2 Materials (in contact with medium)

Sensor element: Platinum and ceramic (zirconium oxide)
Process connection: Stainless steel 1.4435 (AISI 316L)

Only for CLS82E-**NA* 1) and CLS82E-**NB* 2):

Seal: EPDM

- 1) Connection DN25 standard
- 2) Connection DN25 B. Braun

10.5.3 Surface roughness

 $R_a < 0.38 \ \mu m$

Memosens CLS82E Index

Index

Endress+Hauser

A	Process
Ambient temperature	Process pressure
	Process temperature
C	Product identification 6
Charles 16	Product safety 5
Check Connection	R
Mounting	Repair
Connection	Repeatability
Check	Response time
Ensuring the degree of protection 11	Return
D	S
Degree of protection	Safety
Ensuring	Operation
Technical data	Product
Disposal	Workplace safety
Documentation	Safety instructions
Е	Sensor
Electrical connection	Calibration
Environment	Cleaning
I	Connecting
	Spare parts
Incoming acceptance	Storage temperature
Intended use	Surface roughness
M	Т
Materials	Technical data
Maximum measured error 16	Environment
Measured variables	Input
Measuring ranges	Mechanical construction 18
Measuring uncertainty	Performance characteristics 16
Mounting 10	Process
Check	Temperature compensation
N	remperature, pressure raungs
Nameplate 6	U
0	Use
Operational safety	W
Operational safety	Warnings
P	Weight
Performance characteristics 16 Pressure/temperature ratings 18	Workplace safety 4

19



www.addresses.endress.com