# Operating Instructions Raman Rxn2





## **Table of Contents**

T	About this document4	ł
1.1	Warnings	ł
1.2	Symbols on the device	ł
1.3	U.S. export compliance	ł
1.4	Glossary	5
2	Basic safety instruction7	7
2.1	Requirements for the personnel	7
2.2	Designated use	7
2.3	Electrical safety	7
2.4	Operational safety	7
2.5	Lifting provisions for non-cart-mounted Raman Rxn2	3
2.6	Product safety	3
2.7	Important safeguard	3
2.8	Health and safety considerations	)
2.9	Safety and handling notice	)
2.10	Laser safety	)
3	Product description12	2
<b>3</b> 3.1	Product description   12     The Raman Rxn2 analyzer   12	2
<b>3</b> 3.1 3.2	Product description    12      The Raman Rxn2 analyzer    12      Raman RunTime software overview    13	23
<b>3</b> 3.1 3.2 3.3	Product description       12         The Raman Rxn2 analyzer       12         Raman RunTime software overview       13         Product design       14	2 3
<ol> <li>3.1</li> <li>3.2</li> <li>3.3</li> <li>3.4</li> </ol>	Product description       12         The Raman Rxn2 analyzer       12         Raman RunTime software overview       13         Product design       14         Probe Connectors       16	2 3 4 5
<ol> <li>3.1</li> <li>3.2</li> <li>3.3</li> <li>3.4</li> <li>4</li> </ol>	Product description       12         The Raman Rxn2 analyzer       12         Raman RunTime software overview       13         Product design       14         Probe Connectors       16         Incoming product acceptance and identification       17	2 3 4 5 7
<ol> <li>3.1</li> <li>3.2</li> <li>3.3</li> <li>3.4</li> <li>4</li> <li>4.1</li> </ol>	Product description       12         The Raman Rxn2 analyzer       12         Raman RunTime software overview       13         Product design       14         Probe Connectors       16         Incoming product acceptance and identification       17         Incoming acceptance       17	2234577
<ul> <li>3.1</li> <li>3.2</li> <li>3.3</li> <li>3.4</li> <li>4.1</li> <li>4.2</li> </ul>	Product description       12         The Raman Rxn2 analyzer       12         Raman RunTime software overview       13         Product design       14         Probe Connectors       16         Incoming product acceptance and identification       17         Incoming acceptance       17         Scope of delivery       18	<b>2</b> 2 3 4 5 <b>7</b> 3
<ul> <li>3.1</li> <li>3.2</li> <li>3.3</li> <li>3.4</li> <li>4.1</li> <li>4.2</li> <li>5</li> </ul>	Product description       12         The Raman Rxn2 analyzer       12         Raman RunTime software overview       13         Product design       14         Probe Connectors       16         Incoming product acceptance and identification       17         Incoming acceptance       17         Scope of delivery       18         Installation       19	2 2 3 4 5 7 7 3
<ul> <li>3.1</li> <li>3.2</li> <li>3.3</li> <li>3.4</li> <li>4.1</li> <li>4.2</li> <li>5.1</li> </ul>	Product description12The Raman Rxn2 analyzer12Raman RunTime software overview13Product design14Probe Connectors16Incoming product acceptance and identification17Incoming acceptance17Scope of delivery18Installation19Site requirements19	22345773 3
<ul> <li>3.1</li> <li>3.2</li> <li>3.3</li> <li>3.4</li> <li>4.1</li> <li>4.2</li> <li>5.1</li> <li>5.2</li> </ul>	Product description       12         The Raman Rxn2 analyzer       12         Raman RunTime software overview       13         Product design       14         Probe Connectors       16         Incoming product acceptance and identification       17         Incoming acceptance       17         Scope of delivery       18         Installation       19         Site requirements       19         Initial hardware setup       20	<b>2</b> 2 3 4 5 <b>7</b> 7 3 <b>)</b> 9 0
<ul> <li>3.1</li> <li>3.2</li> <li>3.3</li> <li>3.4</li> <li>4.1</li> <li>4.2</li> <li>5.1</li> <li>5.2</li> <li>5.3</li> </ul>	Product description12The Raman Rxn2 analyzer12Raman RunTime software overview13Product design14Probe Connectors16Incoming product acceptance and identification17Incoming acceptance17Scope of delivery18Site requirements19Initial hardware setup20Initial Raman RunTime set up25	22345773995
<ul> <li>3.1</li> <li>3.2</li> <li>3.3</li> <li>3.4</li> <li>4.1</li> <li>4.2</li> <li>5.1</li> <li>5.2</li> <li>5.3</li> <li>6</li> </ul>	Product description12The Raman Rxn2 analyzer12Raman RunTime software overview13Product design14Probe Connectors16Incoming product acceptance and identification17Incoming acceptance17Scope of delivery18Installation19Site requirements19Initial hardware setup20Initial Raman RunTime set up25Electrical and I/O connections28	2 2 3 4 5 7 3 <b>)</b> 9 0 5 <b>3</b>

7	Commissioning	29
7.1	Automation connection	29
7.2	Electrical and I/O connections	31
7.3	Interior of the Raman Rxn2	37
7.4	Raman Rxn2 hardware components	38
8	Operation	40
8.1	Raman RunTime embedded software	40
8.2	Calibrations	41
8.3	Data collection	46
8.4	Using models	53
8.5	Data transfer	56
8.6	Automation connection	58
8.7	Software security	59
9	Diagnostics and troubleshooting	63
9.1	Diagnostics	63
9.2	Warnings and errors	66
9.3	OPC connection troubleshooting	69
9.4	Restarting the system	69
9.5	Raman Rxn2 system and power loss	70
9.6	Recovery console	70
10	Maintenance	73
10.1	Optimization	73
10.2	Servicing the Raman Rxn2 analyzer	75
10.3	Support	77
11	Repair	78
11.1	Servicing and spare parts	
12	Technical data	80
12.1	Specifications	80
12.2	Phazard Area Certification	82
13	Supplementary documentation	83
14	Index	84

## **1** About this document

## 1.1 Warnings

Meaning
This symbol alerts you to a dangerous situation. Failure to avoid the dangerous
situation can result in a fatal or serious injury.
This symbol alerts you to a dangerous situation. Failure to avoid this situation
can result in minor or more serious injuries.
This symbol alerts you to situations which may result in damage to property.

Table 1. Warnings

## 1.2 Symbols on the device

Symbol	Description	
	The Laser Radiation symbol is used to alert the user to the danger of exposure to hazardous visible laser radiation when using the Raman Rxn2 system.	
Â	The High Voltage symbol that alerts people to the presence of electric potential large enough to cause injury or damage. In certain industries, high voltage refers to voltage above a certain threshold. Equipment and conductors that carry high voltage warrant special safety requirements and procedures.	
Intertek	The ETL Listed Mark provides proof of product compliance with North American safety standards. Authorities Having Jurisdiction (AHJ) and code officials across the US and Canada accept the ETL Listed Mark as proof of product compliance to published industry standards.	
X	The WEEE symbol indicates that the product should not be discarded as unsorted waste but must be sent to separate collection facilities for recovery and recycling.	
CE	The CE Marking indicates conformity with health, safety, and environmental protection standards for products sold within the European economic area (EEA).	

Table 2. Symbols

## **1.3 U.S. export compliance**

The policy of Endress+Hauser is strict compliance with U.S. export control laws as detailed in the website of the <u>Bureau of Industry and Security</u> at the U.S. Department of Commerce.

## 1.4 Glossary

Term	Description
AC	Alternating current
ALT	Alternate
ANSI	American National Standards Institute
ATX	Advanced technology eXtended
ATEX	Atmosphere explosible
AWG	American wire gauge
°C	Celsius
CAL	Calibration
CDRH	Center for Devices and Radiological Health
CFR	Code of Federal Regulations
cm	Centimeter
COLL	Collection
CSM	Calibration switching module
CSV	Comma separated value
DC	Direct current
EMC	Electromagnetic compatibility
EO	Electro-optical
EPL	Equipment protection level
EU	European Union
EXC	Excitation
FC	Ferrule connector
GLP	Good laboratory practice
GMP	Good manufacturing practice
HCA	Raman calibration accessory
HPLC	High performance liquid chromatography
Hz	Hertz
IEC	International Electrotechnical Commission
INTLK	Interlock
IP	Internet protocol
IPA	Isopropyl alcohol
IR	Infrared
IS	Intrinsically safe
LED	Light emitting diode
LVS	Low voltage safety
mm	Millimeter
MPE	Maximum permissible exposure
MT	Mechanical transfer

Term	Description
mW	Milliwatt
NA	Numerical aperture
NAT	Network address translation
nm	Nanometer
NOHD	Nominal ocular hazard distance
OPC	Open Platform Communications
PAT	Process analytical technology
РСМ	Power control module
PDF	Portable document format
PLC	Programmable logic controllers
p/n	Part number
QbD	Quality by design
RTU	Remote terminal unit
SOP	Standard operating procedure
SPC	Spectrum
ТСР	Transmission control protocol
UA	Unified architecture
UDP	User datagram protocol
UPS	Uninterruptible power supply
USB	Universal serial bus
USP	SIMCA project file
V	Volt
W	Watt
WEEE	Waste Electrical and Electronic Equipment

Table 3. Glossary

## 2 Basic safety instruction

### 2.1 Requirements for the personnel

- Installation, commissioning, operation, and maintenance of the measuring system may be carried out only by specially trained technical personnel.
- The technical personnel must be authorized by the plant operator to carry out the specified activities.
- The electrical connection may be performed only by an electrical technician.
- The technical personnel must have read and understood these Operating Instructions and must follow the instructions contained therein.
- Faults at the measuring point may only be rectified by authorized trained personnel. Repairs not described in this document must be carried out only directly at the manufacturer's site or by the service organization.

### 2.2 Designated use

The Raman Rxn2 analyzer is designed for use in the following applications:

- Chemical composition measurements of solids, liquids, gases, or turbid media in a laboratory or process development environment. The Raman Rxn2 is particularly suited for use in the following applications:
  - Endpoint monitoring of chemical reactions.
  - Crystallinity monitoring of solid materials.
  - Critical process parameter monitoring and control in upstream cell culture or fermentation bioprocesses.
  - Molecular structure and composition of plant-based proteins, dairy solids, and cell-based foods.
  - Small molecule pharmaceutical polymorph identification and monitoring.

Use of the device for any purpose other than that described, poses a threat to the safety of people and of the entire measuring system and is not permitted.

## 2.3 Electrical safety

As the user, you are responsible for complying with the following safety conditions:

- Installation guidelines.
- Local standards and regulations electromagnetic compatibility.
- The product has been tested for electromagnetic compatibility in accordance with the applicable international standards for industrial applications.
- The electromagnetic compatibility indicated applies only to a product that has been properly connected.

## 2.4 Operational safety

Before commissioning the entire measuring point:

- 1. Verify that all connections are correct.
- 2. Ensure that electrical cables and optical fiber connections are undamaged.
- 3. Do not operate damaged products, and protect them against unintentional operation.
- 4. Label damaged products as defective.

During operation:

- 1. If faults cannot be rectified: products must be taken out of service and protected against unintentional operation.
- 2. Keep the door closed when not carrying out service and maintenance work.

#### 

#### Activities while the analyzer is in operation introduce risk of exposure to measured materials.

- Follow standard procedures for limiting exposure to chemical or biological materials.
- Follow workplace policies on personal protective equipment including wearing protective clothing, goggles and gloves and limiting physical access to analyzer location.
- Clean any spills using the appropriate site policies on cleaning procedures.

#### 

#### Risk of injury from analyzer door stop mechanism

If the enclosure needs to be open, always open the analyzer door fully to ensure the analyzer door stop engages properly.

### 2.5 Lifting provisions for non-cart-mounted Raman Rxn2

The Raman Rxn2 has incorporated a total of four lifting handles – two on each side of the base unit.

#### 

▶ Two people are required to lift and move the Raman Rxn2 instrument.

The Raman Rxn2 is designed to be lifted and moved by two persons in the horizontal orientation. One person must be on each side of the base unit, with each person using both provided handles.



Figure 1. Raman Rxn2 integrated lifting handles

#### 

• Disconnect fiber cables before moving a non-cart mounted instrument.

## 2.6 Product safety

The product is designed to meet state-of-the-art safety requirements, has been tested, and left the factory in a condition in which it is safe to operate. The relevant regulations and international standards have been observed. Devices connected to the analyzer must comply with the applicable safety standards. See Laser Safety  $\rightarrow \square$  for additional details.

## 2.7 Important safeguard

- Do not use the Raman Rxn2 for anything other than its intended use.
- Do not drape the power cord over counters or on hot surfaces.
- Do not open the enclosure of the Raman Rxn2.
- Do not look directly into the laser beam.
- Do not stare or focus a laser in a diffused direction.
- Do not point a laser at a mirrored surface.
- Do not leave attached and unused probes uncapped or unblocked.
- Avoid shiny surfaces and always use a laser beam block.
- Probe shutter on the Rxn-10 probe should always be in the **OFF** position when probe is not in use.

## 2.8 Health and safety considerations

It is the user's responsibility to understand and comply with all applicable safety regulations. These will be variable based on the installation location of the instrument. Endress+Hauser takes no responsibility for determining the safe use of the instrument based on this qualification procedure.

The following actions and laser safety precautions must always be observed while using the Raman Rxn2:

- The Raman Rxn2 is a <u>Center for Devices and Radiological Health</u> (CDRH) Class 3B device. The user should wear appropriate eye protection.
- The Raman Rxn2 should only be used in a location with a suitable and stable power supply.
- If an interlock is required, all entryways to the room or area housing the Raman Rxn2 analyzer must be fitted with warning signs on doors into the Class 3B area.

## 2.9 Safety and handling notice

Raman Rxn2 analyzers incorporate a 532 nm, 785 nm, or 993 nm laser excitation source. Take the following precautions when handling the analyzer and probes when the laser is **ON**:

- Turn **OFF** the laser power (using the key on the front of Raman Rxn2) before making fiber connections and probe inspections.
- If using a Rxn-10 probehead, turn laser output shutter to the **OFF** position before the removal or attachment of optics.
- Do not look directly into the fiber probe output (when the optic is disconnected) or the output (window) of any
  probes. If a Rxn-10 probehead is found with no optic or probe attached Turn laser output shutter on the Rxn-10
  probehead to OFF or the laser power key on the Raman Rxn2 analyzer to OFF.



Figure 2. Rxn-10 probehead showing the ON and OFF shutter positions

#### NOTICE

Handle probes and cables with care. Fiber cables should NOT be kinked and should be routed to maintain minimum bend radii (~6 inches). Permanent damage to the cables may result if they are bent beyond the minimum radius.

## 2.10 Laser safety

Laser light presents special safety hazards not associated with other light sources. All laser users, and others present, need to be aware of the special properties and dangers involved in laser radiation. Familiarity with the Raman Rxn2 and the properties of intense laser radiation will aid in the safe operation of the Raman Rxn2. The Raman Rxn2 may contain an 532 nm, 785 nm, or 993 nm laser. Refer to your system specification information to determine which laser you have. The combination of intense monochromatic light concentrated in a small area means that, under certain conditions, exposure to laser light is potentially hazardous. In workplace environments, a laser safety program provides environmental, training, and safety controls which may reduce the risk of laser-related injuries or workplace damage. For more assistance with taking appropriate precautions and setting the proper controls when dealing with lasers and their hazards, refer to the most current version of <u>ANSI for Safe Use of Lasers Z136.1</u>. The Raman Rxn2 analyzer has hardware safety controls to reduce the risk of laser-based injuries including an interlock and a spring-loaded protective cap covering the laser output of the fiber-optic cables.

The beam is routed from the rear panel of the instrument via a fiber optic patch-cord. In the unlikely event that the fiber optic probe cable is removed, the interlock is overridden, and the spring-loaded protective cap is overridden,

there will be a laser beam exiting the analyzer unit. This beam has a beam diameter of 103  $\mu m$  and a Numerical aperture (NA) of 0.29.

Table 4 provides the fiber core size and mode and the nominal ocular hazard distance equation for the case of laser exiting directly from the analyzer unit.

Base Unit Used	Fiber Core Size and Mode	Nominal ocular hazard distance (NOHD) Equation
Rxn2 Standard	103 µm multi-mode (NA =0.29)	$r_{\rm NOHD} = 1.7/NA \ (\Phi/\pi MPE)^{1/2}$
		multimode equation
MPE at 532 nm continuous viewing: 1 x 10 <sup>-3</sup> W·cm <sup>-2</sup>		
MPE at 785 nm continuous viewi	ing: 1.479 x 10⁻³ W∙cm⁻²	
MPE at 993 nm continuous viewi	ing: 3.854 x 10⁻³ W·cm⁻²	
$\Phi$ = Maximum Power in Watts (V	N)	

Table 4. Laser safety

Another nominal hazard zone calculation needs to be performed to account for the scenario when the analyzer is equipped with a probe. Depending upon the probe utilized, the beam diameter, numerical aperture of the fiber optical cable to the probehead and focusing characteristics of the probehead, the nominal hazard zone calculation will change depending upon if the potential exposure point is at the tip of the probe or at a broken optical fiber. Refer to the specifications section in the pertinent Endress+Hauser Raman probe operating instructions for the appropriate information to complete the nominal hazard zone calculations pertaining to other exposure points.

#### 

Laser beams can cause ignition of certain substances such as volatile chemicals. The two possible mechanisms for ignition are direct heating of the sample to a point causing ignition and the heating of a contaminant (such as dust) to a critical point leading to ignition of the sample.

#### 

- The Raman Rxn2 uses a Class 3B laser as defined in <u>ANSI Z136.1</u>. Direct eye contact with the output beam from the laser will cause severe damage and possible blindness.
- Use of controls or adjustments or performance of procedures other than those specified in this manual may result in hazardous radiation exposure.

For more assistance on appropriate precautions and setting the proper controls when dealing with lasers and their hazards, refer to the most current version of <u>ANSI for Safe Use of Lasers Z136.1</u>.

#### 2.10.1 Optical safety

The Raman Rxn2 is outfitted with a Class 3B laser. Lasers at 785 nm and 993 nm present further safety concerns because the radiation is nearly invisible. Always be aware of the initial direction and possible scattering paths of the laser. The use of OD3 safety glasses is highly recommended for 532 nm and 785 nm excitation wavelengths and OD4 for a 993 nm excitation wavelength.



Figure 3. Laser safety glasses

#### 2.10.2 Electrical safety

The Raman Rxn2 utilizes AC and DC voltages inside the enclosure. Do not disassemble the laser enclosure as there are no serviceable parts inside the laser assembly. Only qualified personnel familiar with high voltage electronics should open the system enclosure to perform necessary maintenance or service.

### 2.10.3 CDRH compliance

The Raman Rxn2 is designed and built to meet the laser performance requirements of 21 <u>Code of Federal</u> Regulations (CFR), Chapter I, Subchapter (J) and is registered with the CDRH.

The product report for the Raman Rxn2 can be found under accession number 1110121-000.

#### 2.10.3.1 Protective housing

The Raman Rxn2 is enclosed in a protective housing to prevent human access in excess of the limits of Class I radiation as specified in U.S. 21 CFR Section 1040.10 (f) (1) except for the output, which is Class 3B.

#### 2.10.3.2 Remote interlock connector

The Raman Rxn2 is supplied with a remote interlock connector for each channel. This connector allows the operator to utilize an external interlock circuit in conjunction with Raman Rxn2 operations. Design and function of an external interlock circuit should meet the capability and intent of the most current revision of the <u>ANSI Z136.1</u> Standard. No laser radiation for a particular channel is emitted unless both the fiber and remote interlock connectors are connected.

#### 2.10.3.3 Key control

The Raman Rxn2 utilizes a key control system. The laser radiation will not be accessible until the system key switch is in the **ON** position. The key cannot be removed when the switch is in the **ON** position.

#### 2.10.3.4 Compliance labels

The Raman Rxn2 analyzer is certified to comply with the U.S. 21 CFR, Chapter I, Subchapter (J), as administered by the CDRH.

#### 2.10.4 WEEE directive compliance

The Raman Rxn2 complies with the <u>Waste electrical and electronic equipment</u> (WEEE) Directive 2012/19/EU. The WEEE Symbol shown below is placed on all WEEE-compliant assemblies.



Figure 4. WEEE symbol

If no other means of disposal are available, Endress+Hauser offers a "Take Back" disposal program at no cost. To participate in the "Take Back" disposal program, refer to our website (<u>https://endress.com/contact</u>) for a list of contacts in your area.

#### 2.10.5 Specific conditions of use

- 1. The fiber optic cable shall be installed so that the minimum bend radius specified by the cable manufacturer is not exceeded.
- 2. Where it is necessary to monitor the process level to ensure that the optical beam is not exposed to a potentially explosive atmosphere, the devices used to monitor the level shall be intrinsically safe or classed as simple apparatus, or be installed so as to provide a fault tolerance of 2 for Equipment protection level (EPL) Ga equipment or a fault tolerance of 1 for EPL Gb equipment. The functional safety of this arrangement has not been assessed as part of this certification and it is the responsibility of the installer/user to ensure that an appropriate mechanism is in place.
- 3. Where Intrinsically safe (IS) galvanic isolators are added to the main enclosure in order to produce IS signals to external apparatus not covered by this certification, the IS galvanic isolators shall have an ambient working temperature upper limit of at least 55 °C (131 °F). The IS parameters pertaining to these isolators shall be conveyed to the user in an appropriate manner. The IS nature of any such circuits has not been assessed as part of this certification and this certificate is not to be taken as indication that these IS circuits comply with relevant requirements.

## 3 Product description

## 3.1 The Raman Rxn2 analyzer

The Raman Rxn2 analyzer, powered by Kaiser Raman technology, is a for-purpose embedded system with built-in Raman RunTime control software. Raman spectroscopy provides the chemical specificity of mid-IR spectroscopy and the sampling simplicity of near-IR spectroscopy. By operating in the visible or near-infrared spectral region, Raman spectroscopy allows vibrational spectra to be collected *in situ*, using fiber-coupled probes, without sample purging, and without the use of speciality sampling devices. There are four possible configurations of the Raman Rxn2 analyzer: single-channel, four-channel, hybrid, and starter. All Raman Rxn2 analyzers employ a unique self-monitoring system to ensure the validity of each analysis. The analyzer is capable of two-point self-calibration in extreme environments and utilizes self-diagnostics and spectral correction methods when system calibration is unnecessary. The analyzer's precision is essential for robust chemometric analyses and calibration transfer between instruments. The Raman Rxn2 suite of analyzers allow for remote fiber-optic connections to probe sampling points for installation flexibility. And, all configurations of the Raman Rxn2 analyzer are designed for use with the Endress+Hauser line of Raman fiber optic probes and optics. An ergonomic mobile cart (including built-in probe and optic storage) is available as an option for all Raman Rxn2 configurations.

### 3.1.1 The Raman Rxn2 single-channel and four-channel analyzer

The Raman Rxn2 single channel analyzer offers one fiber-optic sampling probe connector for the accurate measurement, collection, monitoring, and analysis of a single sample. The Raman Rxn2 four channel analyzer provides four fiber-optic sampling probe connectors. Both the Raman Rxn2 single-channel and four-channel analyzers are designed for use in an analytical or process development laboratory for routine sample measurements, quality assurance, or process development applications in life sciences, chemical and food and beverage industries. The four-channel Raman Rxn2 was designed for customers who need to support process development activities so that they can monitor multiple vessels. The ability to follow several different reactions at the same time can quickly build process knowledge and simplify technology transfer from a laboratory into a process environment. With the Raman RunTime software embedded into the analyzer, the Raman Rxn2 single-channel and four-channel analyzers meet the needs of both Good laboratory practice (GLP) and Good manufacturing practice (GMP) governed areas within the pharmaceutical industry for Process analytical technology (PAT) and Quality by design (QbD) applications.

Raman Rxn2 single channel or four channel analyzers can be configured with a 532 nm, 785 nm, or 993 nm excitation wavelength laser.

### 3.1.2 The Raman Rxn2 hybrid analyzer

The Raman Rxn2 hybrid analyzer is configured with two fiber optic sampling probe connectors: one Rxn-20 probe channel and one Alternate (ALT) non-Rxn-20 probe channel. The Raman Rxn2 hybrid analyzer is a unique Raman analyzer because it features the Rxn-20-based large volumetric probe and a backscattered probe. These two probe types enable a variety of applications for solids, liquids, and turbid media. A backscattered immersion probe is the preferred approach for measuring liquids because of its short focus, optical window, and bubble-shedding design. The Rxn-20 probe is optimized for large volumetric measurements, enabling focus-free, non-contact representative measurements of solids or turbid media. Since the Raman Rxn2 hybrid is equipped with these two probes, this analyzer provides maximal sampling flexibility for laboratory, quality control, and process development purposes. With the Raman RunTime software embedded into the analyzer, the Raman Rxn2 hybrid analyzer meets the needs of both Good laboratory practice (GLP) and Good manufacturing practice (GMP) governed areas within the pharmaceutical industry for Process analytical technology (PAT) and Quality by design (QbD) applications.

The Raman Rxn2 hybrid analyzer is only available with a 785 nm excitation wavelength laser.

### 3.1.3 The Raman Rxn2 Starter analyzer

The Raman Rxn2 Starter analyzer is a complete kit comprised of a lower-resolution single-channel Raman Rxn2 analyzer and a Rxn-10 probehead. While configured for a single probe connector, the Raman Rxn2 Starter can be configured up to four channels. The Raman Rxn2 Starter is designed for cart-portable or benchtop uses such as material quality, reaction monitoring, basic science research, quality assurance, and unknowns identification. The variety of non-contact or immersion optics compatible with the Rxn-10 probehead provides sampling flexibility to support a variety of applications.

The Raman Rxn2 Starter analyzer is only available with a 785 nm excitation wavelength.

## 3.2 Raman RunTime software overview

Raman RunTime embedded software is the control platform for the suite of Raman Rxn analyzers. Raman RunTime software is intended for easy integration with standard multivariate analysis and automation platforms to enable a real-time, *in situ* process monitoring and control solution. Raman RunTime presents an OPC and Modbus interface, which provides clients with analyzer data as well as analyzer control functions. Raman RunTime is fully embedded into Raman Rxn analyzers.

The main view of Raman RunTime displays four quadrants for four channel analyzers, one for each probe and a status bar (bottom) for a quick view of warnings and acquisition status. Single channel analyzers display only one main channel/probe window, while hybrid analyzers display two channel/probe windows, one for the Rxn-20 probe channel and one for the non-Rxn-20 probe ALT. Batch details are accessed and edited from each corresponding quadrant/probe window. To switch back and forth between the main view and batch detail views, click the **Title Bar** for each probe/quadrant. Views of current spectrum vs. process values (model results) can also be easily swapped by clicking on the **Quadrant/Probe** window display. Additional features, such as system settings, calibration and diagnostics, are found under the Options section on the lower left corner of the screen.



Figure 5. Raman RunTime's main view in a Raman Rxn four channel analyzer

#### 

▶ When the Raman Rxn2 analyzer's main power switch and laser key are turned **ON**, probes should be shuttered or covered, or remain immersed in the sample to be measured.

#### NOTICE

• See Operation  $\rightarrow \square$  for more information about working with the Raman RunTime software.

## 3.3 Product design

#### 3.3.1 Front panel

On the front panel of the instrument are the standard user interfaces. These include the main **ON/OFF** power switch, the laser **ON/OFF** key switch, Light emitting diode (LED) indicators, and a Universal serial bus (USB) 3.0 port.



Figure 6. Front panel of a Raman Rxn2 four channel analyzer

#	Name	Description
1	Laser Key Switch	The laser key switch turns the laser on and off. The <b>Red</b> LED indicator adjacent to the laser key switch indicates the laser power status. To activate turn the key to the <b>ON</b> position.
2	Main Power Switch	The main power switch turns the instrument on and off, which includes the laser regardless of the position of the laser key switch. The <b>Power</b> push button incorporates a <b>Blue</b> LED in the shape of a power symbol, which indicates the system power status. The <b>Power</b> push button will communicate error conditions using blink codes when embedded software is not able to communicate them. To turn the instrument on, press and release the <b>Power</b> button once. To turn a responsive instrument off, shut down using Raman RunTime. If the instrument is unresponsive, it may be powered down using a long 10 second press and hold of the <b>Power</b> button.
3	USB 3.0 Port	The USB 3.0 port is intended to obtain diagnostic exports from the instrument using a USB flash drive.
4	Probe Connection Status Indicators	The bank of <b>Yellow</b> LED indicators between the laser key and USB 3.0 port indicate the physical connection status of the probes. While the Raman Rxn2 four channel analyzer front panel has four LED indicators, the front panel of the Raman Rxn2 hybrid analyzer has only two LED indicators, and the front panel of the Raman Rxn2 single channel analyzer has only one LED indicator.

Table 5. Front panel of Raman Rxn2

### 3.3.2 Rear panel

On the rear panel of the instrument are standard Input/output (I/O) ports. These include touchscreen, USB, ethernet, serial, and video ports.



Figure 7. Rear external circuit input/output panel of a Raman Rxn embedded analyzer

#	Name	Description
1	Touchscreen USB Port	USB 3.0 port used to connect to the touchscreen.
2	USB Port (Auxiliary)	USB 3.0 backup port. Reserved for future use.
3	Ethernet Port (2)	Ethernet ports for the network connection.
4	Touchscreen Video Port	Touchscreen video port for connection to local touchscreen display (if needed).
5	RS-485 Serial Port	RS-485 serial port, half-duplex. Provides automation data via Modbus Remote terminal unit (RTU). Port settings configurable in Raman RunTime.

Table 6. Raman Rxn2 ports

#### 3.3.3 Base unit rear panel – single channel and four channel analyzers

All normal system Input/output (I/O) is located at the rear of the base unit. This includes:

- EO fiber connector/electrical connections for up to four remotely installed probes for the four channel Raman Rxn2 analyzer (single channel analyzers have only one probe connection). The electrical connection contained within the fiber optic is an intrinsically safe interlock loop which will turn off the laser in case of fiber breakage.
- Four remote interlock connections for the four channel Raman Rxn2 analyzer (only one for the single channel configuration), each intrinsically safe and in series with the fiber breakage detection loops described in previous bullet.
- Two TCP/IP Ethernet ports for OPC and Modbus automation as well as remote control.
- One serial RS-485 port for Modbus automation.
- One Mini Display Port for local display (if needed).
- Two USB 3.0 Type A ports, one for local touchscreen (if needed) and one reserved for future use.
- AC power inlet, C13 plug required. See Specifications → 
   <sup>(a)</sup>

#### 

Care must be taken to make the bend radii of fiber optic cable breakouts as large as possible (minimum of 6 inches) with NO kinks.

#### 3.3.4 Base unit rear panel – hybrid analyzers

All normal system I/O is located at the rear of Raman Rxn hybrid analyzers. This includes:

- Fiber optic connectors/electrical connection for one remotely installed Rxn-20 probe. The electrical connection contained within the fiber optic is an intrinsically safe interlock loop which will turn off the laser for the Rxn-20 probe in case of fiber breakage. All connections are protected by a cover secured to the rear panel with two socket head cap screws.
- EO fiber connector/electrical connection for one remotely installed ALT probe. The electrical connection contained within the fiber optic is an intrinsically safe interlock loop which will turn off the laser for the alternate probe in case of fiber breakage.
- Two remote interlock connections for the Rxn-20 and ALT probes, each intrinsically safe and in series with the fiber breakage detection loops described in previous bullets.
- Two TCP/IP Ethernet ports for OPC and Modbus automation as well as remote control.
- One serial RS-485 port for Modbus automation.
- One Mini Display Port for local display (if needed).
- Two USB 3.0 Type A ports, one for local touchscreen (if needed) and one reserved for future use.
- AC power inlet, C13 plug required. See Specifications → 
   <sup>(a)</sup>

#### 

• Care must be taken to make the bend radii of fiber optic cable breakouts as large as possible with NO kinks.

### 3.4 Probe Connectors

The probes are connected to the base unit at the connection panel located on the rear of the base unit.

For single channel or four channel Raman Rxn2 configurations, as well as the Alternate (ALT) probe channel of the Raman Rxn2 hybrid configuration, each channel employs a single robust Electro-optical (EO) connector which contains excitation and collection fiber-optics as well as an electrical laser interlock loop. The interlock contained within the probe fiber optic is a low voltage and current loop designed to detect fiber cable breakage and will turn laser emission off for that channel in the event of a break. Ensure latch is engaged after inserting EO fiber connector.

For the Rxn-20 channel of the Raman Rxn2 hybrid instrument, the main fiber optic bundle is furcated into three FC fiber optic connectors and one electrical interlock loop connector. The FC fiber optic connections are used for laser excitation, Raman scatter collection, and auto-calibration. The interlock loop contained within the probe fiber-optic is a low voltage and current loop designed to detect fiber cable breakage and will turn laser emission off for the Rxn-20 probe in the event of breakage.

## 4 Incoming product acceptance and identification

### 4.1 Incoming acceptance

- 1. Verify that the packaging is undamaged. Notify the supplier of any damage to the packaging. Keep the damaged packaging until the issue has been resolved.
- 2. Verify that the contents are undamaged. Notify the supplier of any damage to the delivery contents. Keep the damaged goods until the issue has been resolved.
- 3. Check that the delivery is complete and nothing is missing. Compare the shipping documents with your order.
- 4. Pack the product for storage and transportation in such a way that it is protected against impact and moisture. The original packaging offers the best protection. Make sure to comply with the permitted ambient conditions.

If you have any questions, please contact your supplier or your local sales center.

#### NOTICE

#### Incorrect transportation can damage the analyzer

• Always use a lifting truck or a forklift to transport the analyzer.

#### 4.1.1 Nameplate

The nameplate located on the rear of the analyzer provides the following information about your device:

- Manufacturer Contact Information
- Laser Radiation Notice
- Electric Shock Notice
- Model Number
- Serial Number
- Wavelength
- Maximum Power
- Build Month
- Build Year
- Patent Information
- Certification Information

Compare the information on the nameplate with the order.

#### 4.1.2 Identifying the product

The order code and serial number of your product can be found in the following locations:

- On the nameplate.
- In the delivery papers.

#### 4.1.3 Manufacturer address

Endress+Hauser 371 Parkland Plaza Ann Arbor, MI 48103 USA

## 4.2 Scope of delivery

The scope of delivery comprises:

- Raman Rxn2 analyzer in the configuration ordered
- Raman Rxn2 Operating Instructions
- Raman Rxn2 Certificate of Product Performance
- Local declarations of conformity, if applicable
- Certificates for hazardous zone use, if applicable
- Raman Rxn2 optional accessories, if applicable

If you have any queries, please contact your supplier or local sales center.

## 5 Installation

### 5.1 Site requirements

The base unit enclosure contains all functional components of the analyzer. The enclosure is designed for orientation in either the vertical tower position or the horizontal desktop position. The air flow through the base unit draws air in from the front of the unit and exhausts at the rear of the unit. To limit inlet air obstruction, at least 15.5 cm (6 inches) of clearance must be maintained at the front of the unit. To limit exhaust air obstruction, the rear of the base unit must have 15.5 cm (6 inches) of clearance to the nearest obstruction.

### 5.1.1 Electrical power

The supply voltage is recommended to be regulated and free of voltage spikes. It is recommended, but not required, that an Uninterruptible power supply (UPS) be used with the instrument.

For a Raman Rxn2 analyzer, the location selected must have four (4) voltage power outlets, one (1) for each of the following:

- Base unit.
- Touchscreen monitor.
- Optional power meter or the Raman calibration accessory (HCA).

Alternatively, these components may be plugged into a power strip on one (1) outlet.

#### 5.1.2 Location

The Raman Rxn2 analyzer can be located on a level surface such as a laboratory benchtop or an equipment cart. In addition, the location selected should be:

- Free of moisture, dust, and corrosive vapors.
- Isolated from excess vibrations.
- Protected from direct sunlight.

#### 5.1.3 Ventilation

The location selected should allow for adequate ventilation to both the front and rear of the base unit. A space of no less than 6 inches (152.4 mm) must be provided in both the front and rear of the base unit for proper inlet and outlet air movement.

#### 5.1.4 Temperature

The Raman Rxn2 base unit and touchscreen monitor are designed to operate within a temperature range of 15  $^{\circ}$ C to 30  $^{\circ}$ C (59  $^{\circ}$ F to 86  $^{\circ}$ F). In any installation case, care must be taken to keep the instrument inlet air and surrounding air within this temperature range.

#### 5.1.5 Relative humidity

The Raman Rxn2 base unit and touchscreen monitor are designed to operate within an ambient relative humidity range of 20 % to 80 %, non-condensing.

## 5.2 Initial hardware setup

#### 5.2.1 Installation of Raman Rxn2 analyzer

Endress+Hauser requires in some cases, that installation and initial analyzer setup be performed only by trained service professionals from Endress+Hauser or its affiliated channel partners. The section provides only a basic overview of the analyzer setup process, and does not facilitate a full onsite installation or IQ/OQ. For IQ/OQ, it is required that a trained representative from Endress+Hauser or one of its authorized distributors be present for both initial inspection and installation of the Raman Rxn2 analyzer. Prior to the installation, prepare the site per the requirements in Section 5.1.

#### 5.2.2 Connecting the touchscreen monitor

1. Connect a touchscreen monitor to the **Display** and **Touchscreen USB** ports on the back of the instrument.



Figure 8. Touchscreen display connections

#### 5.2.2.1 For single-channel, four-channel, and Starter analyzers

Raman Rxn2 single-channel, four-channel, Starter, and hybrid (ALT channel only) analyzers utilize a unique fiberoptic connector with integrated electrical laser interlock loop. The EO fiber connector on the Raman Rxn2 analyzer is referred to as the 'device-side' connector. The EO fiber connector on the probe fiber cable is referred to as the 'cableside' connector. The instrument device-side connector uses an integrated spring-loaded cap that will protect the internal fibers from contamination.

To connect a probe:

1. Remove cover from probe fiber cable-side connector.



Figure 9. Electro-optical fiber connection cover

- 2. Clean fiber tips of cable-side connector prior to installation if cleanliness of fiber tips is unknown.
  - First use a lens wipe very lightly saturated with a solvent, such as reagent grade acetone or 100 % Isopropyl alcohol (IPA) followed by a final clean with a 1.25 mm fiber cleaning tool. Do not use the same wipe for both fiber tips.
  - Swipe the fiber tip once with the damp portion of the wipe, then swipe once more with a dry portion of the same wipe. Repeat for both fiber tips.



Figure 10. Cleaning electro-optical fiber connection

3. Next, use an IBC 1.25 mm ferrule cleaner with the bulkhead adapter attached to do a final clean of the center of the ferrule where the fiber resides. Press together until a click is heard and repeat once.



Figure 11. Final clean of electro-optical fiber connector fiber tips

4. Release latch and open spring-loaded cap on the Raman Rxn2 analyzer device-side connector. Insert cable-side connector into instrument device-side connector fully and engage latch to secure. The connectors are polarized and can only be inserted one way. The slotted head screws on the face of both connectors should be facing out.



Figure 12. Connecting an electro-optical fiber cable to a Raman Rxn2 four channel analyzer

5. Repeat for each probe.

#### 

Unused probes that are attached to the Raman Rxn2 analyzer should ALWAYS be capped to prevent stray ambient light from entering the probe. Stray ambient light, whether from an uncapped probe or through incomplete light shielding of the sample, may produce undesirable spectral interferences and result in calibration failure or inaccuracy.

#### 

- Probes attached to the Raman Rxn2 analyzer should always be capped or pointed away from people toward a diffuse target if not installed in a sample chamber.
- 6. For each EO fiber cable, restrain the probe EO fiber optic cable using the strain relief located at the rear of the instrument.

#### 

When installing the probe in situ, the user must provide strain relief to the EO fiber optic cable at the probe installation location.

#### 5.2.2.2 For hybrid analyzers

The ALT channel utilizes an EO cable with an integrated electrical laser interlock loop, and the process of connecting the ALT channel is described in the previous section. The Rxn-20 channel of the Raman Rxn2 hybrid analyzer has four connections points: fiber electrical interlock, excitation, collection, and calibration. The excitation and calibration connections utilize FC style fiber connectors, while the collection connection utilizes a Mechanical transfer (MT) style fiber connector. A remote interlock connector is also provided for the Rxn-20 channel and is located adjacent to the fiber electrical interlock connection. The Rxn-20 excitation, collection, and calibration fibers are fragile, and care must be taken to properly route and restrain the fibers using the following procedures.

To connect a probe to the Rxn-20 channel:

1. Remove the Rxn Rxn-20 plastic cover by unscrewing the two socket head cap screws using the 7/64 in. ball driver included with the instrument. This will reveal the individual connectors for the Rxn-20 channel.



Figure 13. Rxn-20 channel cover on rear of Raman Rxn2 hybrid analyzer and connectors for Rxn-20 channel

- 2. On the cable-side connectors, clean ONLY the excitation and calibration fiber tips prior to installation if cleanliness of these fiber tips is unknown.
  - First use a lens wipe very lightly saturated with a solvent, such as reagent grade acetone or 100 % IPA followed by a final clean with a 2.5 mm fiber cleaning tool. Do not use the same wipe for both fiber tips.
  - Swipe the fiber tip once with the damp portion of the wipe, then swipe once more with a dry portion of the same wipe. Repeat for both fiber tips.
- 3. Employ the two strain reliefs on the rear of Raman Rxn2 hybrid analyzer (using the 7/64 in. ball driver) to secure the fiber cable assembly.



Figure 14. Employing strain reliefs to secure the fiber assembly for the Rxn-20 channel

4. Remove the threaded cap from the Calibration (CAL) port.

5. Connect the clean calibration fiber to the CAL port, aligning the key on the fiber connector with the notch on the CAL port connector. The fiber should be routed in a helical pattern as shown below to avoid kinks.



Figure 15. Proper connection and routing of the calibration fiber

6. Remove the rubber cap from the MT style collection fiber connector. Lift and hold the cover of the Collection (COLL) port as the fiber is connected. Align the white dot on the MT style fiber connector with the white mark on the COLL port and insert the fiber connector until it clicks into place. The fiber should be routed in a helical pattern as shown below to avoid kinks.



Figure 16. Proper connection and routing of the collection fiber

- 7. Remove the excitation fiber clamp by loosening the clamp screw with a 3/32 in. ball driver (provided with analyzer) and sliding it out of the way.
- 8. Remove the threaded cap from the Excitation (EXC) port.



Figure 17. Proper removal of the excitation fiber clamp and threaded cap from the excitation port

9. Connect the clean excitation fiber to the EXC port on the connection panel, aligning the key on the fiber connector with the notch on the EXC port connector. The fiber should be routed in a helical pattern to avoid kinks.



Figure 18. Proper connection and routing of the excitation fiber

- 10. Re-install the excitation fiber safety clamp using a 3/32 in. ball driver.
- 11. Connect the fiber electrical interlock connector to the Interlock (INTLK) port. Route cable behind fiber connections.



Figure 19. Proper connection of interlock connector and fiber routing

12. Replace the Rxn-20 plastic cover, ensuring that no fibers are caught between the cover and connection panel flange. Replace and tighten the two socket head cap screws to hold cover in place using a 7/64 in. ball driver.



Figure 20. Before and after reattaching the Rxn-20 channel c/over

#### 

Unused probes that are attached to the Raman Rxn2 analyzer should ALWAYS be capped to prevent stray ambient light from entering the probe. Stray ambient light, whether from an uncapped probe or through incomplete light shielding of the sample, may produce undesirable spectral interferences and result in calibration failure or inaccuracy.

#### **WARNING**

The beam emerging from the Rxn-20 probe is hazardous to the eye. Always secure the probe so that it is pointing safely away from any personnel. Never handle the probe freely when it is operational.

#### 5.2.3 Powering up the Raman Rxn2

Press and release the **Power** button and turn the laser key switch to the **ON** position. The **Power** button will blink once per second until Raman RunTime launches. The Laser Enable LED will illuminate **Red** and the power switch will illuminate a solid **Blue**.

#### NOTICE

• See Front Panel  $\rightarrow \square$  for more information about the the **Power** button.

#### 5.2.4 Powering down the Raman Rxn2 system

There are two ways to properly power down the Raman Rxn2 analyzer. One of these two methods should always be used to power down the instrument unless it is unresponsive.

1. In Raman RunTime embedded software, navigate to **Options > System > General**. Click **Shut Down**. The analyzer will power down after about 5 seconds.

c	Seneral
INSTRUMENT	
Instrument Name	IOT-Analyzer >
1	Restart
Sh	ut down

Figure 21. Steps to shut down the device via Raman RunTime

OR

1. Press and hold the push button power switch until it starts to blink (2 seconds). Release the push button. A power down message is sent to Raman RunTime, and it will begin power down activities. The instrument will power down after about 5 seconds.

To perform a **hard shutdown**, there are two methods that can be used. These methods should **only** be used if Raman RunTime is unresponsive.

1. Press and hold the push button power switch for at least 12 seconds until the instrument powers down. Then release the button.

#### NOTICE

- After 2 seconds, the power switch will begin to blink; ignore and continue to hold down power button until instrument powers down.
- 2. Unplug the instrument.

NOTICE

▶ See the *Raman RunTime Operating Instructions (BA02180C)* for more information.

### 5.3 Initial Raman RunTime set up

#### 5.3.1 Raman RunTime network configuration

The table below summarizes the ports relevant to specific features. Ports for OPC Classic are assigned dynamically in the indicated range; fixed-port behavior is only available for OPC UA. It is only necessary to open firewall ports for the features intended to use. Ports in bold should be regarded as mandatory for the associated protocol; other ports are recommended for full functionality (e.g. discovery services).

Protocol	Ports
OPC UA	TCP 4840, 4843, 52601, and <b>62886</b>
OPC Classic	TCP <b>135</b> and <b>49152-65535</b>
Modbus	TCP <b>502</b>
File Sharing	TCP <b>139</b> and <b>445</b>
Remote Access	TCP <b>3594</b> and <b>4526</b> for Secure connection; TCP <b>3593</b> , <b>4525</b> , and 5674 for Standard connection; UDP 3702

#### 5.3.2 Initial setup of Raman RunTime

- 1. Customize the analyzer name in **Options > System > General**. For example, Raman Rxn2-785 sn0012345. Click **Apply**. This step is important because the analyzer name is how the system is identified via diagnostic exports.
- 2. Calibrate the touch screen (if needed):
  - Select **Options > System > General > Calibrate Touch Screen**.
  - Follow on screen prompts. A tip for achieving a better calibration is to use the edge of one's fingernail when following on screen prompts and touching the requested touch points.
- Connect to a network using one of the available ethernet ports. View and customize network settings, at Options
   System > Network.
- 4. Set the date and time either manually or via network connection by specifying the time server address at **Options > System > Date & Time**. If setting the date and time manually, first ensure the time zone is set up correctly before proceeding to other adjustments. This step is also important because spectral acquisition and resulting files and OPC tags are managed by the system's date/time.
- 5. Specify names for each probe/quadrant (such as Probe 1) at **Probe Detail view > Settings Tab > Name > Apply**.

#### 

• You should allow the system to stabilize for a minimum of two hours before proceeding to the calibration steps.

#### 5.3.3 Initial calibration and verification

Follow these instructions to calibrate and verify your Raman Rxn2 analyzer for the first time. All calibrations are required prior to spectral acquisition. Raman RunTime will not allow spectra to be collected without passing internal and probe calibrations. Passing the verification step is not required but highly recommended.

Navigate to **Options > Calibration**. The next page shows different calibration window previews for the different Raman Rxn2 analyzer configurations.



Figure 22. Calibration window with quick view buttons for a Raman Rxn2 four channel analyzer



Figure 23. Calibration window with quick view buttons for a Raman Rxn2 single channel analyzer

Calibration		
Internal Calibration	Probe Calibra	tion
Calibrate	1 Calibrate	Verify
2022-06-17 19:38	ON 🚫 2022-06-19 17:15 🚺	not verified
01h	Calibrate	Verify
Calibrate Periodically	UN 💓 2022-00-17 19.32	2022-00-20 10.13
Calibration Report		
View Save		
		Close

Figure 24. Calibration window with quick view buttons for a Raman Rxn2 hybrid analyzer

## 6 Electrical and I/O connections

## 6.1 Overview of rear panel electrical and I/O connections



Figure 25. Rear panel of Raman Rxn2 four-channel analyzer

#	Name	Description
1	Remote Interlock Connectors	Safety feature. To interrupt the laser, remove the black plug.
2	EO Fiber Connector	Provides fiber optic laser radiation output, fiber optic Raman collection and electrical laser interlock loop for each instrument channel. The electrical laser interlock loop is intrinsically safe and is governed by Endress+Hauser drawing 4002396. Match 3 prongs on probe to 3 plugs on EO. Pull latch down to secure probe in place. Laser radiation may NOT exit a channel whose EO fiber connector is removed because removing the EO connector also interrupts the laser interlock loop for that channel.
3	Air Exhaust	Air exhaust fan and outlet.
4	CDRH Product Label	Product information about the Raman Rxn2 analyzer.
5	Analyzer Ports	Touchscreen USB Port, USB Port, Ethernet Ports, RS-485 Serial Port, and Touchscreen Video Port.
6	AC Inlet 100-240 VAC 50/60 Hz	Power socket outlet that AC power to the base unit. Ground pin on this connector serves as the Protective Conductor Terminal.
7	Strain Relief	EO fiber cable strain relief mounting location.

Table 8. Port connection

## 7 Commissioning

### 7.1 Automation connection

Raman RunTime provides network-connected clients with analyzer data as well as analyzer control functions. OPC UA is the recommended protocol because it allows transfer of large data (full spectral data and diagnostics) and is a more reliable connection than OPC Classic. However, legacy support for OPC Classic (DCOM, also called OPC DA) clients is also built-in.

#### NOTICE

The Raman Rxn2 system must be connected to a network for OPC functionality. Network settings can be viewed and configured in **Options > System > Network**.

#### 7.1.1 OPC connectivity

To access the OPC server, all clients must authenticate with the name and password of an authorized user.

- Automation. The built-in user with the name 'kaiser-opc' is always authorized for OPC access. The password for 'kaiser-opc' is 'opc' by default; this password can be changed under Options > System > Automation.
- Security. Additionally, the OPC server can be accessed with user credentials created and managed under Options
   > System > Security.
- OPC UA (recommended). OPC UA clients access the OPC server using UA's binary protocol to connect to
  opc.tcp://<computer-name>, providing the name and password of an authorized user. The "computer name" can
  be viewed or changed under Options > System > Network.

The Basic128Rsa15, Basic256, and Basic256Sha256 algorithm suites are supported for signing and optionally encrypting OPC communications.

 UA Certificates. The first time that a given UA client connects to a Raman RunTime system, it is necessary for that client to accept or "trust" the certificate that the OPC server provides to identify itself. The mechanisms for accepting the server's certificate vary by client.

The OPC server provides a "self-signed" certificate; accordingly, it is common for the client to display a warning that the server's identity cannot be verified. Such a warning is normal, and the server's certificate must be accepted to enable the client to connect.

The OPC server's certificate is tied to the computer name of the Raman RunTime system. Any time the computer name is changed, a new certificate will be generated automatically, and UA clients will need to accept the new certificate before they can connect again.

- OPC Classic (deprecated). Full OPC functionality is available to OPC Classic clients. However, OPC Classic entails
  additional set-up requirements, and many opportunities exist for client-side configuration to interfere with
  successful communication, often in ways that are challenging to diagnose. Also, the underlying technology of
  OPC Classic can limit installation flexibility. For these reasons, it is recommended to use OPC UA instead of OPC
  Classic whenever possible.
- The OPC Classic client program must be executed using credentials that exactly match (name and password) an authorized user on the Raman RunTime system.
- This implies that a Windows user account must be created on the client workstation with a name matching that
  of a Raman RunTime user, and that their passwords must be kept in sync.

#### 7.1.1.1 Asynchronous OPC updates

In addition to responding to polling requests from clients, the OPC server is capable of providing asynchronous updates in which the server notifies clients of changes in the values of subscribed tags without the need for clients to continually poll the server.

#### NOTICE

• Many clients will fail to connect unless asynchronous updates are enabled.

To enable asynchronous updates, the 'kaiser-opc' user must exist on the client workstation with a password that matches the Raman RunTime system, regardless of whether the client program accesses the server using the 'kaiser-opc' user credentials.

#### NOTICE

- ► If the 'kaiser-opc' password is changed under **Options** > **System** > **Automation**, the Raman RunTime system must be restarted under **Options** > **System** > **General** for the new password to take effect.
- Client workstations must reside in the same TCP/IP subnet as the analyzer, with no intervening Network address translation (NAT).

#### 7.1.2 Modbus connectivity

The Raman Rxn2 provides automation data via Modbus TCP (Ethernet) or Modbus RTU (RS-485 serial). If models are running on the instrument, the prediction values available over Modbus may be configured at **Options > System > Automation > Modbus**. In addition to prediction values, other instrument diagnostic values are available over Modbus.

#### NOTICE

See *Raman RunTime Automation Modbus (p/n 4005298)* to view the current Modbus Map for Raman RunTime.

## 7.2 Electrical and I/O connections

An I/O panel is located at the center of the rear panel, which provides several extra low voltage, non-intrinsically safe external circuits:

- Touchscreen. Type A USB 3.0 for connection to a local touchscreen display. This is not required for the
  instrument to operate as all operation parameters may be set through the instrument's automation interfaces. A
  touchscreen kit, which includes the interface cable for this connection, is available for purchase from
  Endress+Hauser under *p/n 2018086*.
- Display. Mini DisplayPort for video connection to a local touchscreen display. This is not required for the
  instrument to operate as all operation parameters may be set through the instrument's automation interfaces.
  This port does NOT support DP++; therefore, an active adapter is needed if connecting to a display without native
  Display Port. A touchscreen kit, which includes the interface cable for this connection, is available for purchase
  from Endress+Hauser under *p/n 2018086*.
- Aux. Type A USB 3.0 reserved for future use.
- **RS-485.** DB9 which provides a half-duplex serial RS-485, two-wire plus ground, Modbus RTU automation interface. Pin 2 is Data +, Pin 3 is Data -, Pin 5 is Ground. All other pins are inactive.

Recommended wiring is commercially available shielded, 2 twisted pair, 22 American wire gauge (AWG), terminated with a DB9 receptacle and back shell kit. Endress+Hauser recommends Carol C1352A cable, TE Connectivity 5-747905-2 receptacle and 1991253-9 back shell kit. Cable and connector/back shell of equivalent specifications may be substituted. One pair is used for Data+ and Data- and one of the wires from the second pair is used for ground. It is not recommended to use the shield as a signal ground. No provision for connecting the shield drain is provided at the Raman Rxn2. The shield may be connected to earth at the device on the opposite end of the cable attached to the Raman Rxn2.

- Network 1. 10/100/1000 RJ45 Ethernet interface. Provides remote control option and automation data via OPC UA, OPC Classic and Modbus TCP. Use standard Ethernet cabling.
- Network 2. Same as Network 1. Both interfaces may be used at the same time.



Figure 26. Rear panel on a Raman Rxn2 four channel analyzer



Figure 27. Rear panel of Raman Rxn2 hybrid analyzer with Rxn-20 channel housing attached (left) and removed (right)

#### 7.2.1 Power and grounding

The Raman Rxn2 has a standard IEC-320 C-14 inlet for power at the rear of the instrument. Any power cord with an IEC-320 C-13 plug will connect to the base unit. The Raman Rxn2 accepts AC power from 100 – 240V and 50/60 Hz. For U.S. applications, a power cord is supplied. For non-U.S. applications, the user must supply a power cable that meets local/national standards.

Also included on the rear of the instrument is a functional earth terminal for additional grounding if required. Primary grounding occurs through the IEC power inlet plug ground terminal which should be connected to the building grounding system.

Do not position the Raman Rxn2 in a way that makes it difficult to remove the mains cord. Only use adequately rated power cables with the Raman Rxn2 system.



## 7.2.2 Electrical interconnect block diagram – single channel Raman Rxn2

Figure 28. Single channel Raman Rxn2



## 7.2.3 Electrical interconnect block diagram – four channel Raman Rxn2

Figure 29. Four channel Raman Rxn2



## 7.2.4 Electrical interconnect block diagram – Raman Rxn2 hybrid

Figure 30. Raman Rxn2 hybrid



### 7.2.5 Hazardous area installation drawing

Figure 31. Hazard area installation drawing

#### 7.2.6 Remote laser interlock connectors

Adjacent to each instrument probe connector resides a remote interlock connector. This connector allows integrators to access the analyzer's individual channel interlock circuit, and connect external switching devices, e.g. emergency stop button, door switch, etc., as additional means to stop laser radiation for each channel. This connector is in series with the associated channel's probe connector interlock loop. The output of this connector is intrinsically safe. Connection of external devices, including cabling, to this connector are governed by drawing 4002396.

The instrument ships with shorting plugs installed in each channel's remote interlock connector. Should field wiring be required for an external switch, a remote interlock pigtail may be purchased to facilitate connection to field wiring using p/n 2018748 (single pigtail) or p/n 2018749 (four pigtails). Replacement shorting plugs may be purchased using p/n 2018429.

If one emergency switch is required to interrupt all four channels of a four-channel instrument simultaneously, a 4pole switch shall be used. The four separate interlock loops are electrically isolated and may NOT be tied together electrically. Endress+Hauser recommends IDEC XN1E-BV404MR for a 4PST-NC emergency stop button switch.

Refer to drawing 3000095 for remote interlock connection details.
## 7.3 Interior of the Raman Rxn2

The interior of the Raman Rxn2 with the cover removed is shown below. The internal components are common among all configurations.



Figure 33. Embedded controller locked in service position

## 7.4 Raman Rxn2 hardware components

### 7.4.1 Laser

The laser used in the Raman Rxn2 is a special version of Endress+Hauser's Class 3B laser. It has a lockable injector that requires no routine adjustment.

The laser can be turned on and off using the LASER ENABLE switch on the front of the base unit at any time while the Raman Rxn2 is operating.

#### 7.4.1.1 Laser apertures

Apertures for laser emission are found in the following places on and within the Raman Rxn2:

- Probehead
- Connection Panel
- Calibration Module
- Laser

#### 7.4.1.2 Laser interlock circuit

If a fiber breaks the interlock, the laser output for that channel is disabled. Other channels with intact fiber interlocks will continue to have the laser available to them.

For laser output to occur on a channel, a probe and remote interlock shorting plug, Endress+Hauser p/n 2018429, must be installed to their respective connectors.

Auxiliary interlock laser emission indicators are on the probes. See the specific probe operating instructions for more information.

## 7.4.2 Spectrograph

The spectrograph includes the optical elements used to filter out the Rayleigh scatter and focus the Raman scatter onto the detector. The spectrograph assembly (which includes the detector) in the Raman Rxn2 is contained in a sealed assembly and has no user-serviceable parts.

## 7.4.3 Calibration switching module

The Calibration switching module (CSM) is a key component of the Raman Rxn2. Using high-reliability switches, it routes the several channels and performs automatic wavelength calibration with a neon source and laser wavelength calibration with an internal Raman shift standard. It also contains a software-controllable shutter for the laser.

The fluorescent lamp provides a broad spectrum of lines for wavelength calibration across the Raman spectrum. The neon lamp also is not susceptible to spectral shift with changes in temperature or pressure as are calibration protocols that rely on Raman bands.

The calibration can be performed on the internal neon lamp without reconfiguring the analyzer to accommodate an external calibration unit. The internal Raman shift standard allows the laser wavelength to be tracked.

Because the calibration light source is internal to the Raman Rxn2, it is susceptible to interference from stray light entering the connected probes. Prevent stray light from entering any probe connected to the base unit by covering the ends of the attached probes.

#### NOTICE

► See Calibrations  $\rightarrow \triangleq$  for more information on calibrating the Raman Rxn2. Intensity calibration requires Endress+Hauser's *Raman Calibration Accessory (p/n 2004909-501)*.

### 7.4.4 Fuses

There are no fuses to replace on the Raman Rxn2. The Raman Rxn2 is powered by a special adapted Advanced technology eXtended (ATX) power supply, which has no external fuses. If a short circuit were to occur inside the Raman Rxn2 it would be on the DC output side of the power supply. In this event, the power supply would turn itself off and would require a manual power reset (i.e., removing the plug for five (5) minutes) after the source of the short circuit was resolved.

### 7.4.5 Mobile cart accessory option

An ergonomic mobile cart/trolley is available as an option for all Raman Rxn2 configurations. The Raman Rxn2 with mobile cart model includes the Raman Rxn2 embedded analyzer, built-in probe and optic storage, a routine-analysis sample compartment, and fiber storage. For connecting main power to the mobile cart, an integrated internationally approved pigtailed power cord is provided.

The pigtail incorporates a standard IEC-320 C-14 inlet to which any standard and locally approved power cord with an IEC-320 C-13 plug may be attached to provide mains power to the cart accessory. The mobile cart accepts AC power from 100 – 240V and 50/60 Hz. Internal to the mobile cart, all mains power is distributed to the Raman Rxn2 base unit, touchscreen monitor and optional Calibration Accessory using an internationally approved power distribution strip with internationally approved cordage.

For mobile cart options sold within the continental United States, an approved 3.0-meter power cord is provided with the Mobile Cart for attachment to mains power. For mobile cart options sold outside the continental United States, a power cord is NOT included. It is the responsibility of the end user or local distributor to provide a locally approved power cord for attachment to mains power.

#### NOTICE

▶ See the Raman Rxn New Cart Operating Instructions (BA02175C) for more information.

### 7.4.6 Air filter

The Raman Rxn2 incorporates a tacked polyester spun air filter element to reduce dust intake into the base unit. The air filter is accessed by a magnetically secured access panel on the front of the instrument. The air filter should be cleaned with compressed air once every month or if the embedded software is reporting an internal over-temperature error (if ambient temperature is within specification). In extremely dusty conditions, the air filter should be cleaned more often. The air filter has a **Blue** tacky side which should be oriented toward the outside of the base unit.

If a replacement air filter is needed, refer to our website (<u>https://endress.com/contact</u>) for the list of local sales channels in your area.



Figure 34. Filter access

## 8 Operation

## 8.1 Raman RunTime embedded software

Raman RunTime is the embedded control software installed on all the Raman Rxn2 analyzers. It is intended for easy integration with standard multivariate analysis and automation platforms to enable a real-time, *in situ* process monitoring and control solution. Raman RunTime presents an OPC and Modbus interface which provides clients with analyzer data as well as analyzer control functions.

Depending on the instrument configuration being used, the main view of Raman RunTime displays four "quadrants" or windows (one for each of the four probes for the four channel analyzer), two windows (one for each of the two probes of the hybrid analyzer), or one window (for the single probe used in the single channel analyzer). The main view of Raman RunTime also displays a status bar (bottom) for a quick view of warnings and acquisition status. Details are accessed and edited from each corresponding quadrant.



Figure 35. Raman RunTime main screen for a Raman Rxn2 four channel analyzer

To switch back and forth between the main view and detail views, simply click the **Title Bar** for each probe/window as applicable. Views of current spectrum vs. process values (model results) can also be easily swapped with a click on the display windows. Additional features, such as system settings, calibration and diagnostics, are found under the Options section. Raman RunTime is also used as the interface between the Raman Rxn2 analyzer and third-party PAT software suites such as synTQ and SiPAT.

#### NOTICE

This section provides details on configuring and using the Raman Rxn2 with Raman RunTime, but it is not intended to replace the full *Raman RunTime Operating Instructions (BA02180C)* or analyzer installation and training from Endress+Hauser representatives.

## 8.2 Calibrations

Raman Rxn2 analyzers have internal calibration standards for spectrograph and laser wavelength. This section assumes the instrument is being setup for the first time and utilizes the Recalibrate All option, which is only required at initial start-up or when the instrument is moved or serviced.

There are several internal calibration options to choose from:

- Automatic. If the instrument is already calibrated, this will compare the current analyzer response to calibration specifications and will recalibrate if the spectrograph wavelength or laser wavelength is out of spec. If the analyzer is uncalibrated, this will perform an alignment calibration, followed by a full wavelength calibration and full laser wavelength calibration.
- **Recalibrate X Axis.** Forces full wavelength and laser calibrations without first checking whether the analyzer is within spec.
- **Recalibrate All.** This will cause the alignment calibration to be repeated prior to performing full spectrograph wavelength and laser wavelength calibrations. Note that when "Recalibrate All" completes, the intensity calibrations and verifications of all probes will be invalidated.

Steps for Internal Calibration:

- 1. Ensure all probes are capped.
- 2. Click **Calibrate** under Internal Calibration.

For initial startup, click **Recalibrate All** from the Calibration Mode drop-down list and click **Calibrate**.

#### NOTICE

• The **Recalibrate All** option is only required at initial start-up or when the analyzer is moved or serviced.

Calibrate Instrument			]
Calibration Mode			
Recalibrate All		÷	
	Calibrate	Cancel	

Figure 36. Recalibrate all

- For subsequent calibrations, select **Automatic**.
- 3. Select **Calibrate Periodically** and set the frequency.
  - Setting is configurable in one-hour increments.
  - At least once per day (1d 00h) is recommended.

	Internal Calibration
	Calibrate
0	2019-03-01 09:38
_	
	1d 00h
✓ C	alibrate Periodically

Figure 37. Periodic calibration setting

A0048761

## 8.2.1 Probe calibration – single channel, four channel, or hybrid analyzers

A Calibration accessory (HCA) with an appropriate optic adapter is required for probe calibration of single channel, four channel, and hybrid (ALT channel only) Raman Rxn2 analyzers. For the Raman optic system for single use, a single use calibration adapter is also needed.

Ensure the HCA calibration has been calibrated within 1 year and has an accompanying Intensity Reference source spectral file (\*.spc).

Probe calibrations can be executed during active experiments, e.g. if a probe needs to be set up while another probe probe is active. When a probe calibration is triggered, any acquisitions currently in progress will be aborted automatically, and the calibration will proceed. Upon completion of the calibration, active probes will automatically resume normal operation.

Steps for probe calibration:

1. Connect the HCA to the first (or only) probe. Use the set screw to hold the probe in place.



Figure 38. Example of the Rxn-10 probe with a bIO Optic connected to the HCA

- 2. Turn on the HCA and set the switch to Intensity. The intensity indicator on the HCA will blink until the bulb warm up time is complete (15 minutes). Do not proceed until indicator stops blinking.
- 3. For Rxn-10 probes, be sure the laser shutter is set to **ON** (indicator parallel to probe body).
- 4. Loading the HCA Intensity Reference spectral file (\*.spc):
  - Attach the USB memory stick that came with the HCA kit to the USB Extension or USB port on the front of the analyzer. This is only required for new device setup or when the HCA is serviced or changed.
  - For single channel or hybrid (ALT channel) analyzers, click Load on the Probe Calibration section of the screen to browse to the correct Intensity Reference spectral file (\*.spc) located on the memory stick. Click Select File.
  - For a four channel analyzer, click **Calibrate** for the targeted probe, which will result in a pop-up window that lets the user load the Intensity Reference spectral file as above.

Probe Calibration for Channel 1
INTENSITY CALIBRATION
<ol> <li>Present the Calibration Accessory to the probe.</li> <li>Activate 'INTENSITY' on the Calibration Accessory.</li> <li>Click 'Calibrate'.</li> </ol>
Intensity Reference
SN0070001 2019-05-28 · Load Save
Cancel Calibrate

Figure 39. Intensity calibration window in Raman RunTime and memory stick provided with the HCA

- 5. Click **Calibrate** to start the calibration process on the selected probe:
  - **Four Channel Raman Rxn2 System.** Make sure the **Blue** indicator is active on the probe intended for calibration cancel and recalibrate if needed. A countdown timer will start.



Figure 40. Four channel calibration

• **Single Channel Raman Rxn2 System.** There is no **Blue** indicator and instead only a countdown timer will start, indicating probe calibration has begun.



Figure 41. Single channel calibration

 Raman Rxn2 Hybrid System. Make sure the Blue indicator is on the ALT probe. Cancel and recalibrate if needed. A countdown timer will start.



Figure 42. Hybrid calibration

6. Repeat the probe calibration for each active probe (if applicable).

NOTICE

► For analyzers with any unused channels/probes, the warning indicator will turn **Yellow** indicating there are uncalibrated channels. To resolve this erroneous warning, turn off unused probes/channels.

## 8.2.2 Probe calibration – hybrid (Rxn-20 channel) analyzers

For the Rxn-20 channel on a Raman Rxn2 hybrid analyzer, follow the steps below:

- 1. Connect the appropriate optic (3mm, 6mm) to the Rxn-20 probe.
- 2. Connect the appropriate extension tube to the optic. The 3mm optic uses the 3 in. long extension tube, while the 6mm optic uses the 8 in. long extension tube.
- 3. Connect the HCA lamp head to the extension tube. Use the thumb screw to hold the HCA lamp head in place.



Figure 43. HCA lamp head and extension tube attached to Rxn-20 non-contact optic and probe

4. After powering on the HCA, set the switch to Intensity. Let the white light bulb warm up for approximately 15 minutes.

#### 

For HCAs delivered after approximately 01-Nov-2017, the HCA intensity indicator light will blink until the halogen bulb warm-up period is complete. Wait to proceed with probe calibration until the light stops blinking.

- 5. Loading the HCA Intensity Reference spectral file (\*.spc):
  - Attach the USB memory stick that came with the HCA kit to the USB port on the front the analyzer. This is only required for new device setup or when the HCA is serviced or changed.
  - On the probe calibration section of the screen, click **Load** to browse to the correct Intensity Reference spectral file (\*.spc) located on the memory stick. Click **Select File**.

	Calibration
Internal Calibration Calibrate 2020-08-27 19:46 01h Calibrate Periodically	Probe Calibration  PhAT  Calibrate  Aug-27 19:53  Not verify  ALT  Calibrate  Verify  Aug-31 15:10  Not verified
Calibration Report View Save	Intensity Reference: HCA-SN0000115 Calibration Date: 2013-01-17 Load Save
	Close

Figure 44. Calibration window for a Raman Rxn hybrid analyzer

#### 6. Click Calibrate.

Make sure the **Blue** indicator is on the Rxn-20 probe. Cancel and recalibrate if needed. A countdown timer will start.

NOTICE

• For analyzers with any unused channels/probes, the warning indicator will turn **Yellow** indicating there are uncalibrated channels. To resolve this erroneous warning, turn off unused probes/channels.

### 8.2.3 Probe verification

For any Raman Rxn2 analyzers (single channel, four channel, starter, or hybrid), probe verification will verify the calibration results using a standard reference sample, such as 70 % IPA or cyclohexane. This step is not required to collect a Raman spectrum, but it is highly recommended.

#### 

Do not use cyclohexane for verification with any bIO-Optic, Rxn-45 and single-use adapters. Bioprocessing probes such as bIO-Optic, Rxn-45 and single-use adapters, should ONLY use 70 % IPA. Only 70 percent by volume (%v/v) will work. Endress+Hauser recommends using CiDehol 70 by Decon Laboratories. All other probes can use cyclohexane.

For bioprocess probes, use a bIO Sample Chamber for verification:

- 1. Connect the bIO sample chamber at the white Teflon end to the first probe and fill it about half-way with 70 % IPA.
- 2. While holding the attached sample chamber for support (to keep it upright), tilt the probe/optic to a 15-30° angle and tap the sample chamber to ensure any bubbles in the 70 % IPA rise to the top.



Figure 45. Sample chamber

- 3. Click **Verify** under the Probe Calibration section and select 70 % IPA as the verification standard.
- 4. Repeat the probe verification for each active probe.

For all other (non-bioprocess) probes use a sample container appropriate for the probe connection:

- 1. Click Verify under the Probe Calibration section and select cyclohexane or 70 % IPA as the verification standard.
- 2. Repeat the probe verification for each active probe.

How to interact with the probe verification results:

- 1. Access the **Spectrum Viewer** button to display the current verification spectra and allow for addition of previous spectra.
- 2. Click Spectrum Viewer to view the current verification spectrum.

2022-06-19 09:55:35	Analyze	Verificati	on Channel 4
Selected Spectr	a		1
2022-06-19 09:55:35	$\Theta$		
2022-06-17 14:34:20	Θ		
	*		
Exp	Add		1.
n l n	11		
500 1000	1500 2000	2500	3000

Figure 46. Probe verification spectrum viewer

- 3. Click the spectrum's name at the top of the window (click the **Date** and **Time** of the spectrum which appears in **Blue**). Click **Add** to include previous spectra on the plot.
- 4. Spectra can also be exported as \*.spc files (to a USB memory stick) from this view.
- 5. Additionally, the symbols under the **Calibrate** and **Verify** buttons can be clicked for details about the recent calibration/verification.



Figure 47. Probe calibration details

6. Click **Export** to save the verification spectrum to an external USB memory stick that can be inserted in auxiliary USB ports on the front and rear of the device. A network connected drive can also be used for \*.spc file export.

## 8.2.4 Calibration and verification reports

1. On the lower left of the screen under **Calibration Report**, click **View** to see the calibration and verification report. A summary of the internal calibration report appears at the top of the screen.

Cal	ibration ]	Report	Endress+Hauser 🖾
Instrument Name	Raman Analy	zer	
Model	Rxn 785		
Serial Number	123456789		
Channel 1: Stream Channel 2: Stream Channel 3: Stream Channel 4: Stream	A B C D		
Channel	Intensity	Verification	
1	PASSED 2022-06-20 12:21	N/A	
2	N/A	N/A	
3	N/A	N/A	
4	N/A	N/A	
In	ternal Calibratio	DB	
Alignment	20	PASSED 122-06-20 12:18	
Wavelength	1	PASSED	
			Close

Figure 48. Calibration and verification report

- 2. The full calibration and verification report can be viewed on screen by clicking the **Black Arrow** above the **Close** button to scroll down.
- 3. Click **Close** at the bottom of the report window to exit.
- 4. Click **Save** to save the full calibration and verification report locally as a \*.pdf file (to a USB memory stick). The calibration report can also be accessed remotely via network Raman RunTime connection or via OPC.

## 8.3 Data collection

There are several collection modalities available in Raman RunTime including Manual, Continuous, Periodic, Focus, and Snapshot.

Exposure Length and Count (sometimes referred to as accumulations) need to be specified for each collection option, except Focus. Focus may be used to determine exposure length.

- **Exposure Length.** How long the laser will be exposed to the sample.
- **Count.** The number of times the exposure will be repeated for a single spectrum.

The goal of determining an appropriate length and count combination is to maximize the signal-to-noise ratio within the time allotted for sampling without overexposing the detector.

- Ideal detector saturation is between 40 % 70 %.
- Avoid detector saturation below 10 % and above 80 %.

### 8.3.1 Notes about data collection with Raman RunTime

- Batch settings for Collection Mode and Exposure Settings are remembered. This makes it very easy to start a new batch that is like the previous.
- It is not necessary to use the same collection mode for each probe. Collection mode differences and how they
  affect sampling sequence are further described in corresponding sections below.
- Active batches can be paused or stopped at any time using the corresponding buttons.
  - Pausing a batch can be used to not disturb the sample/spectrum while interrupting the experiment, such as exposing a reactor to room light or extracting a sample.
  - $\circ$  ~ It is best to pause at the start of a new sample since the current spectrum/sample will terminate.
  - When collection is paused or stopped, Raman RunTime will automatically proceed to the next probe in sequence (or wait until the next periodic acquisition for another probe). Stopping acquisition of a Raman RunTime batch means the same experiment cannot be restarted new spectra will need to be stored under a new "batch" name and collection of spectra. If there is a chance the user might need to return to the experiment, be sure to pause rather than stop the acquisition process.
- Exposure settings can also be changed during an active experiment. It is not necessary to pause to change
  exposure settings. New settings will take effect on the next sample acquisition.

#### 8.3.2 Dark exposures

Dark exposures are used to reduce noise in spectra by accounting for the noise arising from the detector. For example, 532 nm and 785 nm detectors naturally have low noise, so dark exposure requirements for these systems are minimal and mostly automatic. However, 993 nm detectors have intrinsically high noise levels and require extra dark exposure attention.

Dark handling for Raman Rxn2 532 nm and 785 nm analyzers:

 There is only one user initiated dark setting for 532 nm and 785 nm systems, located by clicking Advanced under Exposure Settings.

Exposure 9	Set	tings		
10.0s	×	75	Advanced	
length		count		A00487

Figure 49. Exposure settings

Leave the Force New Dark check box deselected which should be the default setting for 532 nm and 785 nm systems.



Figure 50. Force new dark

Selecting the **Force New Dark** check box will collect and apply a new dark every time a spectrum is collected. Therefore, this is not recommended for most applications since it will double the collection interval time.

- Dark exposures are automatically collected at the beginning of acquisition whenever an Exposure or Count setting has been changed.
  - There is an exception if a dark exposure has already been collected and stored for the same Length and Count setting, e.g. for another probe or during a previous experiment. Then the system will use the stored dark.
  - To expire an old dark and force a new one at any length and count interval, select the **Force New Dark** check box for at least one interval, then turn it off.
- If a dark exposure is needed when acquiring in continuous or periodic mode, one will be collected at the start of the first acquisition and used for all subsequent spectra.
  - This will cause the first spectral collection interval to take double the time of subsequent intervals.
  - The status indicator specifies when a dark exposure is active and how long is remaining.



Figure 51. Dark exposure

#### NOTICE

It is best to collect a dark after the system has stabilized for about 2 hours after start up. To collect a dark to be stored and used repeatedly, such as for routine exposure settings, allow the analyzer and laser to remain on for 2 hours prior to dark collection if the laser has been turned off. If the analyzer has been relocated, then it should stablize for at least one day with the laser on before collecting a new dark. Note that this is not applicable to moving short distances on a Raman Rxn2 cart.

Dark handling for Raman Rxn2-993 nm single channel and four channel analyzers:

- On 993 nm systems, dark exposures are collected automatically at the end of every acquisition.
- The number of dark exposures collected at the end of each acquisition is determined by the Dark Exposures setting under the **Advanced** button.

- Typically, dark subtraction is effective at mitigating the non-sample contributions in the measurement due to "dark current" when the Dark Exposures setting is approximately one half the count setting.
- In processing the final spectrum, the "trailing" dark exposures taken after the sample is measured are combined with "leading" dark exposures taken before the sample is measured.
- Typically, the trailing dark exposures from one acquisition are used as the leading dark exposures for the next
  acquisition. This significantly reduces the time needed to collect the requisite number of dark exposures for a
  given acquisition.
- Alternatively, new leading dark exposures are collected at the beginning of every acquisition if any of the following conditions are met:
  - $\circ$   $\;$  The length setting has changed since the most recent dark collection.
  - $\circ$  The most recent dark collection completed more than 10 minutes prior.
  - $\circ$  ~ The Force New Dark option is selected under the Advanced button.
- Due to the conditional nature of leading dark collection, there is the potential for inconsistent acquisition times, especially on the first acquisition after starting a new experiment or rebooting the instrument, or generally when using periodic or manual collection. Selecting the Force New Dark option can be used to achieve consistent acquisition times, when doing so is more beneficial than minimizing acquisition time.

## 8.3.3 Focus mode

Focus mode is used to position probes (typically non-contact optics) or determine appropriate collection settings for a given sample or process. No spectra are saved from Focus mode – focus spectra are for visual evaluation only.

#### NOTICE

- ► For the Rxn-20 probe with the Raman Rxn hybrid analyzer, be sure to use the appropriate extension tube for the Rxn-20 optic to help in locating the sample to the focal plane.
- 1. Position probe to sample and click **Focus** under the probe details view.

A Suggested Exposure will be displayed. The example below shows 1.7s – 3.0s for a sample of 70 % IPA.



Figure 52. Focus mode example

2. Move the **Detail** bar to the "slow" position to increase the focus time as needed. Samples yielding weaker signal features may not be obvious with the "fast" focus setting.

The current signal vs. max signal is displayed so that the current probe/sample positioning can be compared to the max observed signal. The current signal will change with repositioning of the probe/sample.

- 3. Reposition the probe or sample to achieve the best focus/signal-to-noise.
- 4. Click **Reset** to reset the remembered max signal.
- 5. Click **Stop** to exit focus mode.

## 8.3.4 Snapshot mode

Snapshot mode is similar to Focus, in that no spectra are saved. Snapshot is intended for quick acquisitions to measure laser power, try out different length and count settings for a sample, or more commonly to quickly assess spectral quality either visually or by applying a model.



Figure 53. Snapshot mode and options

- 1. To access Snapshot mode, click anywhere within the spectral plot from one of the probe detail views. The spectra overlay plot/Snapshot view will appear.
- 2. Select an exposure length and count, then click **Acquire**:
  - The snapshot spectrum will be overlaid with the most recent spectrum if an experiment is already in progress, such as shown below.
  - It will not be stored with the current batch data; a snapshot spectrum is for onscreen tests only.
- 3. Spectra from previous batches/experiments can be overlaid with the snapshot spectrum by clicking the spectrum name (the Blue date and time appearing in the top right header). Then click **Add** and browse to select the reference spectra:
  - Multiple spectra can be added to the overlay, but only a single spectrum is distinguished as the "principal" spectrum by displaying it in **Blue** and listing its name in the header.
  - When a new snapshot is collected, it automatically becomes principal, so is shown in blue. The spectrum is principal and can be changed by selecting/highlighting it in the spectrum list.
  - This list is viewable by clicking **Primary File Name** (located at the top, to the left of **Analyze** button) to display the Selected Spectra drop-down list.
- 4. Additionally, a model can be applied by clicking **Analyze** and selecting a model. Ad hoc analysis only applies to the principal spectrum.



Figure 54. Analyze function in snapshot mode

It is recommended to use Snapshot mode for *in situ* verification of starting materials, especially for long experiments like bioprocesses. Using this method, verification spectra can be collected quickly without needing to adjust batch exposure settings. To exit the spectra overlay view, click anywhere on the quadrant name or click **Exit** in the upper left of the display window.



Figure 55. Exit button to exit the spectra overlay

## 8.3.5 Manual mode

Use the Manual collection mode to acquire spectra that are triggered manually by clicking **Acquire**. A batch/experiment name can be specified to group manual samples together for spectra storage and sample names can be specified for each spectrum.

1. First create a new name by clicking **Edit > New** from the Probe Detail view – the example screen shown at right shows the manual batch name as "Off-line Samples."



Figure 56. Manuacollect mode options

- 2. Set the Collection mode to **Manual**.
- 3. Click **Play**. The Acquisition Batch will show "In Progress" but no spectra will be collected yet until triggered (by clicking **Acquire** in a later step below).
- 4. Use the **Up** or **Down** arrows to adjust Exposure Settings set the Exposure Settings, both length and manual collect mode options count.



Figure 57. Exposure settings as described with length and count boxes

5. If a custom sample name is desired, click the + option next to **Acquire** to enter a sample name, such as the Sample 1 example shown here:

Enter Mar	nual /	Acquisitio	n Name							
Cancel	Sa	mple 1							Acquire	
q	v	v e	e	r	t y	y	u	i	o p	
a	1	s	d	f	g	h	j	k	1	
123		z	x	с	v	b	n	m	back	2
ABC					space				shift	

Figure 58. Keyboard that pops up to rename the sample

- 6. Click **Acquire** to collect a spectrum.
  - To collect another spectrum, specify a new sample name (if desired), change Exposure Settings as needed, then click **Acquire** again.
  - Or click the **Stop** symbol to end the sequence of manual samples stored together.

### 8.3.6 Continuous mode

- Continuous collection mode cycles through each active probe as quickly as possible. It is recommended for method development when collecting frequent reference samples, and for monitoring and controlling when models are active.
- Continuous mode includes a buffered acquisition option.
- Buffered acquisition provides a way to receive updated spectra and model values more frequently, without sacrificing accuracy or repeatability.
- The buffered acquisitions setting specifies the number of consecutive acquisition cycles to combine in each reported spectrum by adding together the individual acquisition spectra.
- The spectrum and model values are updated at the end of each acquisition cycle.
- Only the combined spectrum is reported, the individual acquisition spectra are not.
- Active models operate on the combined spectrum, not on the individual acquisitions.
- Setting buffered acquisitions to a value of 1 (the default) effectively disables the feature, resulting in conventional acquisition in which each spectrum is independent of previous spectra.
- To configure buffered acquisition that achieves the same total target collection time as non-buffered acquisition, reduce the count setting and increase the buffered acquisitions setting, keeping their product the same. For example:
  - If exposure settings of 10s x 60 with 1 buffered acquisition are used, spectra and process values will update only at the end of the full spectrum collection cycle, approximately every 10 minutes.
  - But if settings of 10s x 6 with 10 buffered acquisitions are used, then spectra and process values will update approximately every 1 minute.
  - Each reported spectrum will still include 10 minutes of total acquisition time, because each update adds together the ten most recent acquisitions of 10s x 6.
- For applications requiring long collection intervals, buffered acquisition is not recommended when more than one probe is active.
  - As collection alternates among multiple probes, combining consecutive acquisitions on a given probe runs the risk of including outdated process spectra if other active probes have long intervals.
- 1. To set up a continuous batch, first create a batch under **Edit > New** and specify a new batch name.
- 2. Set the collection mode to **Continuous**.
- 3. Set the exposure and buffer settings desired.



Figure 59. Collection mode details for continuous mode

10068003

- 4. Repeat for all active probes and click **Play** in the desired probe sequence.
  - Pausing, stopping, adding new batches, rebooting, etc. could affect the probe sampling sequence in continuous mode. If cycle order matters, periodic mode is recommended.

## 8.3.7 Periodic mode

Periodic collection mode acquires spectra at specified intervals and is typically used during method development to allow spectra to sync with samples being drawn from reactors or other timed process events. Buffered acquisitions are not an option in periodic mode.

- In the example shown here, periodic acquisition is scheduled for 60 min.
- The batch details show when the next acquisition is due and how many counts (spectra) have been collected for the current batch.
- If spectral acquisition is active in periodic mode, the Next Acquisition Due will read "now."
- 1. To set up a periodic batch, first create a batch under **Edit > New** and specify a new name.
- 2. Set the collection mode to **Periodic**.

Collection Mod	e
Continuous	5
	buffered acquisitions
Periodic	60m 00s
	acquisition period
• Manual	🖸 Acquire   +
Exposure Setti	ngs
10.0s × 7	75 Advanced
length co	ount

Figure 60. Periodic mode batch details

- 3. Set the exposure settings and acquisition period.
- 4. Repeat for all active probes and click **Play** in the desired probe cycle order (1, 2, 3, 4 or 1, 3, 2, 4, etc.).

## 8.4 Using models

Raman RunTime supports executing models generated in SIMCA<sup>®</sup> (<u>Sartorius</u>), GRAMS IQ<sup>™</sup> (<u>Thermo Fisher</u> <u>Scientific</u>), PEAXACT (<u>S-Pact</u>), Unscrambler<sup>™</sup> (<u>Camo Analytics</u>, <u>An Aspen Technologies Company</u>), Solo (<u>Eigenvector</u>), and Data Library. These multivariate models are typically based on in-line Raman spectra correlated to off-line analytical measurements and can be used to turn raw spectra into meaningful process values in real time. Model results can be viewed on-screen as well as communicated via OPC and Modbus.

Raman RunTime supports the following model file versions:

- SIMCA<sup>®</sup> version 13, 14, 15, 16, and 17
- GRAMS IQ<sup>™</sup> versions 9.3 and previous
- PEAXACT version 4, 5, 5.4, 5.6, and 5.7
- Unscrambler<sup>™</sup> version 11
- Solo/PLS\_Toolbox version 8.9 and 9.0

### 8.4.1 Recommendation for model and process value names in Raman RunTime

- Model titles and Y variable names ("Process Values" in Raman RunTime) should be changed in the final models to reflect the desired display names and OPC tags.
- It is wise to include units (g/L, mM, etc.) in the Process Variable name, since model outputs/Raman RunTime do not automatically display units.
- For OPC communication it is very important to use a consistent naming scheme, because each time a Model Title
  or Assay name is changed, the OPC tags will change accordingly (and need to be reconfigured for OPC
  communication).
- The Data Library can be used to ensure consistent Process Value naming. Contact Endress+Hauser for more information on Data Library.
- Avoid using periods and commas in Model Titles and Assay names.
- Special characters (e.g. &, %) work with some OPC configurations, but should always be tested and removed from Model/Process Value names as needed.
- There can be multiple models per Process Value type. This allows for easy prediction comparisons during model development/validation phases, and the use of multiple Process Value models per batch, for example Gluc M1, Gluc M2 or VCD Day 1-6, VCD Day 7-14.

## 8.4.2 Loading models

- 1. Save the desired project files from the modeling package to a USB memory stick and connect it to the USB port on the front of the Raman Rxn2. Alternately, models can be loaded from a network connected folder.
- 2. Select **Options > System > Analysis > Add Model** and browse to the appropriate model file. Repeat for additional files.
- 3. The file name(s) will appear in the Analysis window. Select a file name to turn model components and properties on/off. Repeat for additional files.
- 4. Return to the main view and go to the details section for one probe.
- 5. Press the **Analysis** tab in the lower left and enable the model file(s) on or off. Repeat for additional probes.



Figure 61. Analysis screen to select models

## 8.4.3 Viewing model results

1. From the main view during active experiments, process values will automatically be displayed instead of the current spectra.



Figure 62. View of spectrum for probe 1

- 2. To view the most recent spectrum instead of process values from the main view, click the quadrant display.
  - For hybrid analyzers, the current spectrum and process values are both automatically displayed, there is no need to click the window to change views.
  - For single channel analyzers, there is no quadrant view. Process values will be displayed under the Analysis tab.
  - Model results will update each time a new spectrum is complete, or after each buffer cycle in continuous mode.
  - If models are not active for a probe/batch, then "No Analysis" will appear when the spectral plot is clicked.



Figure 63. Results for a Rxn2 four channel analyzer with models active for probe 1 and probe 4

A0049241 A0049242

To access additional model result information (which is the only way to view model results for single channel analyzers):

- 1. From a probe details view, click the **Analysis** tab.
- 2. Use the **Expand** and **Collapse** buttons to show more or less model details.



Figure 64. Probe details view of model process values and properties

- 3. Click **Trends** to enter the trend viewer.
- 4. Select a **Principal Trend** and, if desired, a **Comparison Trend**.

			Stream	С		
	Principal Trend:		Batches:		Comparison Trend:	
$(\times)$	Bio HT - Glucose (g/L)		✓ Stream C Batch	~	Bio HT - Lactate (g/L)	Ý
15						
10						
5						
0						
-5						
	14:30	14:40	14:50	15:00	15:10	15:2
	Jun 17, 2022					
	Options 🕨					
	man Analuzor		ОК			3 3 4
ГКа	man Analyzer				2022-06-20	10:35:27

Figure 65. Model results in the trend view

### 8.4.4 Saving model results

Stored model results are included in exported batch data and are also available via OPC if Client connections are established.

## 8.5 Data transfer

Raman RunTime offers two ways to transfer or export spectral batch data in \*.spc format. One method is manually and locally exporting on to a USB memory stick. The second is automatically (once configured) exporting \*.spc files over the local network.

## 8.5.1 Batch data export

Spectra (\*.spc format) and model results (\*.csv format) for collected batches can be exported locally from the Raman Rxn2 onto a USB Memory Stick.

- 1. Connect a USB memory stick to the USB port on the front of the Raman Rxn2.
- 2. From the probe detail view, click **Edit**, check the batches to export, then click **Copy**.
- 3. Browse to the USB memory stick and desired folder, and click **Select Folder**.

#### NOTICE

► Two \*.csv files (processvalues.csv and allproperties.csv) containing model results will also be included in the export, ONLY if models were active and predicting during collection.

#### 8.5.1.1 SPC file network export

If the Raman Rxn2 is connected to a network, Spectrum (SPC) files are automatically published in network-accessible folders. A common file share is available on the Raman Rxn2 at \\<*computer-name>\DataLibraryBatchExport*. This directory contains distinct subfolders corresponding to each acquisition batch:

- 1. Batch folder names are of the form *Probe* <#>\_<*batch*>\_<*YYYYmmdd-HHmmss*>, where <#> is the instrument collection channel, <*batch*> is a user-specified batch name, and <*YYYYmmdd-HHmmss*> encodes the date and time the batch was started.
- 2. Spectrum file names are of the form *<batch>\_YYYYmmdd-HHmmss>*, which encodes the time the individual acquisition completed.

## 8.5.2 Remote access

Remote Access permits the Raman RunTime user interface to be accessed using a desktop web browser from a remote workstation, that meets the following requirements:

- Google Chrome and Microsoft Edge are supported web browsers.
- Network connectivity between the instrument and the remote workstation.

There are two ways to access the Raman RunTime user interface remotely:

- **Standard.** Navigate to *http://<IPaddress>:3593* or *http://<hostname>:3593* in a qualifying web browser to use Standard remote access.
- Secure. Navigate to *https://<hostname>:3594* in a qualifying web browser to use Secure remote access. The client must be able to resolve the analyzer hostname for secure connections. Raman RunTime employs a dynamically generated self-signed TLS server certificate in support of encryption. Clients must accept and trust this certificate to successfully communicate with Raman RunTime.

The IP Address and hostname of the analyzer can be viewed and configured under **Options > System > Network**.

When first connecting through secure remote access, clients may be presented with a privacy warning. This warning is expected because the analyzer's self-signed certificate is not trusted by default. The remote client can bypass the warning to access the analyzer by following the steps shared below.

To eliminate the certificate warning from a particular remote workstation, trust can be granted to the analyzer by adding its certificate to the "Trusted Root Certification Authorities" store on the remote workstation. Instructions for this follow.

#### NOTICE

- When Security within Raman RunTime is enabled, an attempt to access Standard remote access will automatically be redirected to the Secure remote access URL.
- A new TLS server certificate will be generated each time the analyzer's hostname is changed.

### 8.5.3 Status indicator

The circular **Status** button in the center of the main screen has three modes: **Green**, **Yellow**, and **Red**. When the status indicator is **Green** with an OK in the center, it means there are no detectable problems.



Figure 66. Green OK status indicator

When the status indicator turns **Yellow** with a flashing warning in the center, it means there is a caution. Click the **Yellow** status indicator for an explanation for the warning.



Figure 67. Explanation for why system has issued a yellow warning

#### NOTICE

► For analyzers with any unused channels/probes, the warning indicator will turn Yellow indicating there are uncalibrated channels. To resolve this erroneous warning, unused channels can be turned off.

When the status indicator turns **Red** with an Error in the center, it signifies a more serious problem that needs immediate attention. Click the **Red** status indicator for an explanation of the error. The error message will disappear automatically once the problem is resolved.



Figure 68. Explanation for why system has issued a red error

Endress+Hauser

## 8.6 Automation connection

Raman RunTime provides network-connected clients with analyzer data as well as analyzer control functions. OPC UA is the recommended protocol because it allows transfer of large data (full spectral data and diagnostics in this case) and is a more reliable connection than OPC Classic. However, legacy support for OPC Classic (DCOM, also called OPC DA) clients is also built-in.

#### NOTICE

The system must be connected to a network for OPC functionality. Network settings can be viewed and configured in **Options > System > Network**.

## 8.6.1 Basic OPC connectivity

To access the OPC server, all clients must authenticate with the name and password of an authorized user.

The built-in user with the name 'kaiser-opc' is always authorized for OPC access. The password for 'kaiser-opc' is 'opc' by default; this password can be changed under **Options > System > Automation**.

The OPC server can be accessed with user credentials created and managed under **Options > System > Security**.

- OPC UA (Recommended). OPC UA clients access the OPC server using UA's binary protocol to connect to opc.tcp://<computer-name>, providing the name and password of an authorized user. The "computer name" can be viewed or changed under Options > System > Network.
- **UA Certificates.** The first time that a given UA client connects to a Raman Rxn2 system, it is necessary for that client to accept or "trust" the certificate that the OPC server provides to identify itself. The mechanisms for accepting the server's certificate vary by client.

The OPC server provides a "self-signed" certificate; accordingly, it is common for the client to display a warning that the server's identity cannot be verified. Such a warning is normal, and the server's certificate must be accepted to enable the client to connect.

The OPC server's certificate is tied to the computer name of the Raman RunTime system. Any time the computer name is changed, a new certificate will be generated automatically, and UA clients will need to accept the new certificate before they can connect again.

OPC Classic (Deprecated). Full OPC functionality is available to OPC Classic clients. However, OPC Classic entails
additional set-up requirements, and many opportunities exist for client-side configuration to interfere with
successful communication, often in ways that are challenging to diagnose. Also, the underlying technology of
OPC Classic can limit installation flexibility. For these reasons, it is recommended to use OPC UA instead of OPC
Classic whenever possible.

The OPC Classic client program must be executed using credentials that exactly match (name and password) an authorized user on the Raman Rxn2 system. This implies that a Windows user account must be created on the client workstation with a name matching a Raman Rxn2 user, and that their passwords must be kept in sync.

## 8.6.2 Asynchronous OPC updates

In addition to responding to polling requests from clients, the OPC server is capable of providing asynchronous updates in which the server notifies clients of changes in the values of subscribed tags without the need for clients to continually poll the server. Note that many clients will fail to connect unless asynchronous updates are enabled.

To enable asynchronous updates, the 'kaiser-opc' user must exist on the client workstation with a password that matches the Raman Rxn2 system, regardless of whether the client program uses the 'kaiser-opc' user credentials.

#### NOTICE

- ► If the 'kaiser-opc' password is changed under **Options > System > Automation**, the Raman Rxn2 system must be restarted under **Options > System > General** for the new password to take effect.
- Client workstations must reside in the same TCP/IP subnet as the instrument, with no intervening NAT.

## 8.6.3 Modbus

The Raman Rxn2 provides automation data via Modbus TCP (Ethernet) or Modbus RTU (RS-485 serial). If models are running on the instrument, the prediction values may be configured at **Options > System > Automation > Modbus**. In addition to prediction values, other instrument diagnostic values are available over Modbus.

## 8.7 Software security

Security can be enabled to block non-authorized users from performing Raman RunTime functions. By default, security is set to **OFF**.

- 1. To turn security settings on browse to **Options > System > Security**.
- 2. Move the Security setting to **ON** and select **OK** for the Administrator password notice.
- 3. Enter a password for the Administrator user and then verify the password in the second password step. Be sure to record the Administrator password. It is required to turn on/off security and create new users.



Figure 69. Setting up security

4. Once the Administrator password is verified, the system security will switch to **ON**.

#### NOTICE

- The Administrator password can be provided to all users, or multiple users with different access levels can be added.
- 5. Click **New** to add a new user.
- 6. Create and verify the new user's password.
- 7. Specify the **Privilege Level** for the user by selecting **User**, **Operator**, or **Administrator** from the drop-down list. See below for privilege level definitions. Once security is **ON**, access to restricted functionality will require the name and password of an authorized user.
- 8. To lock the system immediately after enabling security, click **Lock System Now**.
- 9. Click **OK** when the Lock System? confirmation message appears, and then **OK** again when the System locked message appears.

Upon returning to the main Raman RunTime quadrant/probe window view, a **Blue** lock symbol will appear over restricted features. On this main Raman RunTime view, note that the system can also be locked by clicking **Lock** in the lower left corner next to the Raman analyzer name.

#### NOTICE

When the security feature within Raman RunTime is enabled, the system will automatically lock itself after 15 minutes of inactivity.



Figure 70. Security settings, adding users, and restricted access

## 8.7.1 User privilege levels

The Raman Rxn2 analyzer with Raman RunTime presents a customized embedded spectroscopic platform that is not compatible with installation of third-party software or customer reconfiguration beyond that which is available through the Raman RunTime user interface.

The customized embedded nature of the Raman Rxn2 presents a non-standard attack surface. Raman RunTime's forpurpose operational profile eliminates risky user behaviors and network accessibility is limited to instrument core services. The Raman Rxn2 employs password protection (detailed below) of instrument functionality and only authenticated updates digitally signed by Endress+Hauser can be applied. Furthermore, there is no facility for running non-Endress software, and any system modifications are automatically purged upon restart.

Security can be enabled to block non-authorized users from performing Raman RunTime functions. Security is set to **OFF** by default. There are three user privilege levels: **User**, **Operator**, and **Admin**. The User level is the most basic, followed by Operator with more privileges, and then Admin which can perform all actions. These privileges are detailed below:

User Level	Action
User	<ul> <li>Change subject (channel) display name</li> </ul>
	<ul> <li>View summary and detail of active batch acquisitions</li> </ul>
	<ul> <li>View calibration of information</li> </ul>
	<ul> <li>View diagnostics</li> </ul>
	<ul> <li>Perform exports (Basic, Diagnostic, Full)</li> </ul>
Operator	User level actions plus:
	<ul> <li>View active batch acquisitions</li> </ul>
	<ul> <li>Perform calibration and view calibration information</li> </ul>
	<ul> <li>Perform verification and view verification results</li> </ul>
	<ul> <li>Start/stop batch acquisitions</li> </ul>
	Focus
	<ul> <li>Snapshot</li> </ul>
	<ul> <li>Ad hoc analysis</li> </ul>
	Change system display name
	Configure date and time
	Restart
Admin	No Restrictions
	Operator level restrictions plus:
	<ul> <li>Add/remove models</li> </ul>
	<ul> <li>Configure models (display name, enable/disable components and properties)</li> </ul>
	Configure network
	<ul> <li>Configure Open platform communications (OPC)</li> </ul>
	<ul> <li>Change security settings and manage users</li> </ul>
	<ul> <li>Apply embedded software updates</li> </ul>

Table 9. User privileges

## 8.7.2 Password management and issues

It is a common best practice to routinely change user passwords. In order to do so, an assigned user with varying privilege levels can click **Change Password** in the bottom left corner of the login keyboard screen. Raman RunTime system administrators can also change their passwords in the same way.



Figure 71. Enter user password screen and enter system administrator password screen

A0049216 A0049217

To recover a lost or forgotten password for an assigned user, a user will have to contact their Raman RunTime system administrator to help reset their password. The system administrator will be able to do so by navigating to **Options > System > Security**, then by logging into their system administrator account on this screen. From here, they can select one of the individual assigned users and select **Change Password** at the bottom of the screen. Note than from this screen's Change Password request, it will not ask for the user's old password, just a new password.

Furthermore, if the Raman RunTime system administrator forgets their own password, they can select the **I Forgot** icon in the top right corner of their system administrator password entry screen.

After doing so, they will be provided with a code and will be instructed to contact Endress+Hauser Support and share this code. In response, they will receive a reset code to enter at their prompted screen. They can then select **Authenticate** and proceed to set their new password on the screen that follows:

Reset Password for System Administrator				
ENTER RESET CODE				0
Share the following code with Endress+Hauser Support:	А	В	с	D
GCAHHD	E	F	G	н
Then, enter the reset code provided above.	I	J	к	back
			Cancel A	uthenticate

Figure 72. Password recovery screen for a Raman RunTime system administrator

## 8.7.3 Additional security features

In addition to creating users for password protection of system functionality, the Raman RunTime system has additional security features to protect against unauthorized usage, such as:

- Customized embedded system presents non-standard attack surface.
- For-purpose operational profile eliminates risky user behaviors.
- Network accessibility limited to analyzer core services.
- Only authenticated updates digitally signed by Endress+Hauser can be applied.
- No facility for running non-Endress software.

Unauthorized system modifications are automatically purged upon reboot.

## 8.7.4 Export

An Export file is necessary to restore the instrument, if necessary. A file can be saved to the instrument, networked remote computer or to a flash drive. Go to **Options > Diagnostics > Export**.



Figure 73. Steps to export a file

If the user chooses to create an export file by clicking **New**, they have a choice between **Basic**, **Diagnostic**, and **Full**. Click **Export** to create the file.



Figure 74. Menu received to export new file

The **Basic** button contains the configuration and calibration information, while the **Diagnostic** button provides logs to assist in troubleshooting with Endress+Hauser. The **Full** button contains all comprehensive acquisition and analysis information in addition to information available in Basic and Diagnostic sections.

If chosen, the **Save** option either lets you save the created export file to a USB flash drive or download it to the network computer accessing the embedded RunTime software.



Do you want to open or save Raman Analyzer (20190924-0913).zip from 10.139.70.61? Open Save 
Cancel ×
Figure 75. Steps to save export file

## 8.7.5 Client system requirements

- Google Chrome or Microsoft Edge (for remote access)
- OPC UA client software (for OPC interface)
- OPC Classic client software (for OPC interface)
- Modbus client software (for Modbus interface)

# 9 Diagnostics and troubleshooting

## 9.1 Diagnostics

System diagnostics can be viewed from **Options > Diagnostics > Environment** tab. The purpose of the system's Diagnostics section is for the user to view system environment data, trend historical environment data, and to save an export file to perform a system restore, if necessary. Regular system exports should be done to ensure the system configurations and calibrations are backed up for future use.

NOTICE

- System exports are intended for Endress+Hauser Support to investigate problems or restore analyzers. The contents are subject to change with new RunTime versions and should not be treated as user accessible files.
- Full exports can be very large. A Full export is recommended for regular archival of system info and logs and is required for a system restore.





Figure 76. Steps to reach environment tab and export tab

## 9.1.1 Trends

1. To view system diagnostics, go to **Options > Diagnostics** from the main Raman RunTime menu. This will bring you to the Environment tab within the Systems Diagnostics dialog box where you can see system data.

Name	Value		
stem Environment			
CSM Diamond Heater Temperature	63.0		
Detector Temperature	-40.0		
External Temperature	22.9		
Internal Temperature	24.3		
Laser Cooler Current	0.4		
Laser Cooler Voltage	0.9		
Laser Diode Current	0.9		
Laser Diode Temperature(C)	19.0		
Laser Diode Voltage	2.4		
Laser Enclosure Temperature(C)	22.8		
Laser Heat Sink Temperature(C)	22.9		
Laser Power	333.0		

Figure 77. Environment tab to view system data

2. To display Trends of diagnostic values, click the **Trends** tab. Within the Trends tab, users can access the Principal Trend, History, or Comparison Trend as shown below.



Figure 78. Choices with the trends tab



Figure 79. More choices with the trends tab

## 9.1.2 Diagnostic exports

It is an Endress+Hauser advised best practice to export user/spectral data, custom settings and configurations frequently. A full system Export is also recommended to be generated and saved onto external media periodically. This should be considered for inclusion in the local site Standard operating procedure (SOP).

#### NOTICE

A Full level Export file is required to fully restore the system (including settings, configuration, models, and data) to an earlier state, if necessary.

On the system interface, go to **Options > Diagnostics > Export** tab. To create a new export file, click **New**. The user can opt between Basic, Diagnostic, and Full exports. Click **Export** to create the file.



Figure 80. Steps to create a new export file

#### NOTICE

While the last generated export file is saved onto the onboard analyzer storage, it is only saved there until a newer export file is generated which overwrites the old one. As such, the generated export files should be periodically backed up onto a USB drive for backup and restore purposes.

The Basic option contains the configuration and calibration information, while the Diagnostic option provides logs to assist in troubleshooting with Endress+Hauser. The Full option contains all comprehensive acquisition and analysis information in addition to information available in the Basic and Diagnostic options.

🕦 Stream A 🗖	Stream B <a></a>	System Diagnostics	
Net Started	Dargent 01:42:07	Raman Analyzer (20200922-1928).rex	
3 Stream C	I Stream D		
no analysis	no analysis		
Diagnostics System Options Raman Analyzer	Net Started	Environment Trends Export Close	A0049218

Figure 81. Steps to save an export file

After creating a new export, it will be saved and held on the onboard analyzer storage until overwritten by a newly created one. The Save option on the same screen allows you to either save the created export file to a USB drive or download it to the network computer accessing the remote interface of the Raman RunTime software.

To perform a diagnostic export:

- 1. Navigate to the **Options > Diagnostics > Export** tab. Click **New**.
- 2. Insert a USB flash drive into the USB port on the front of the Raman Rxn2 analyzer.
- 3. Select **Basic**, **Diagnostic**, or **Full** (descriptions are shown on screen).

Full exports are usually very large in file size. A Full export is recommended for regular archival of system information and logs. A Full export may be requested when working with Endress+Hauser Support.

4. Once the export archive has been generated, click **Save** to save the created export file to a USB drive or download it to the network computer accessing the remote interface of the Raman RunTime software.

## 9.2 Warnings and errors

When the system is fully calibrated and operating as expected, the **Status** button in the middle of the main view status bar will read OK and appear **Green**.



Figure 82. Status bar

Symbol	Description
Warning	If a system warning is encountered, the display will change to <b>Yellow</b> . Warnings should be acknowledged but immediate action may not be necessary. Click the <b>Status</b> button to view details of the warning. The most common warning occurs when all the channels are not occupied. The button will pulse continuously until the problem is resolved. Click the <b>Status</b> button to view details about the warning.
Error	If a system error is encountered, the display will change to <b>Red</b> . An error requires immediate action to restore system performance. Click the <b>Status</b> button to view details about the error.

Table 10. Warning and errors

In some cases, users may choose not to utilize all available channels on a Raman Rxn2 analyzer. These unused/uncalibrated channels may result in warnings generated, thereby putting the whole system in a warning state (yellow indicator). To resolve these erroneous warnings about unused channels not being calibrated, the user can individually turn off unused probes/channels in the **Options > Calibration** screen and selecting the **ON/OFF** marker beneath each probe's number as shown below.





	Calibration	1 Stream A 🗖	💿 Stream B 🗖
Internal Calibration Calibrate 2020-09-23 16:41	Probe Calibration 1 Calibrate Verify	no analysis	no analysis
01h Calibrate Periodically	2 Calibrate Verify ↓ ON Ø Sep-23 20:55 ▲ Sep-23 21:01	Net Started 00:00:00 Accession from	Net Started
	Channel Disabled	Channel Disabled	Channel Disabled
Calibration Report View Save	4 Channel Disabled	Options 🕨	
	Close	Raman Analyzer	A0 2020-09-23 21:13:26

Figure 84. Channels 3 and 4 shown disabled and warning indicator resolved to green OK state

If a system error is encountered, the display will change to **Red**.

1. Click the **Red** status indicator to view details about the warning or error.



Figure 85. Analyzer status details

2. In the event the analyzer stops communication with the interface, go to **Options**, select **System**, choose **Restart** and the analyzer will reboot. This will re-establish camera/interface communication.

## 9.2.1 Low laser power

To check laser environment data, go to the **Options > Diagnostics > Environment** tab.

If it is suspected that laser power is low due to lower signal in spectra, conduct a power check with the laser power meter. A laser should register between 380-400 mW of Laser Power. The Laser Diode Current should be 0.7 to 2.0A and will creep up over time due to normal diode ageing.

The user will get a warning from RunTime once laser diode current exceeds 1.6A recommending service of laser module soon, to avoid downtime. Once the laser diode current reaches the 2.0A limit, the laser is considered to be in a failure state, and as a result laser power will gradually start decreasing until it gets to zero. For Technical Service, refer to our website (<u>https://endress.com/contact</u>) for the list of local sales channels in your area.

System Diagnostics				
Name	Value			
System Environment				
CSM Diamond Heater Temperature	64.9			
Detector Temperature	-40.0			
External Temperature	23.0			
Internal Temperature	30.0			
Laser Cooler Current	1.1			
Laser Cooler Voltage	1.6			
Laser Diode Current	0.8			
Laser Diode Temperature(C)	18.9			
Laser Diode Voltage	1.7			
Laser Enclosure Temperature(C)	29.9			
Laser Heat Sink Temperature(C)	31.8			
Laser Power	397.0			
Environment Trends	Export			
		Close		

Figure 86. System diagnostics environment tab

## 9.2.2 ON/OFF button flashing

The **ON/OFF** button emits a flashing blink code to communicate a problem when the software is unavailable.

Sign	Problem	Solution
2 blinks in quick succession followed by a solid long pause.	Indicates problem with main power supply. Could indicate power has been interrupted. Flashing will cease when reserve power is depleted if it is not replenished.	Check security of power cord and its connection. If there is no facility power outage, problem could be with power unit and a replacement is necessary, call service.
6 blinks in quick succession	It is too hot inside the instrument. The device is rated for an environment up to $35 ^{\circ}C (95 ^{\circ}F)$ . Instrument will shut down its power source when it is too hot.	Check ambient temperature in the environment around the instrument. If temperature does not exceed limit, call service.

9.2.3	RunTime system	warnings and erro	rs
-------	----------------	-------------------	----

Diagnostic Title	Faults when
Internal Calibration	Internal calibration missing on multi-channel instrument.
Wavelength Calibration	Wavelength calibration missing on single-channel instrument.
Laser Calibration	Laser calibration missing on single-channel instrument.
Intensity Calibration	Intensity calibration missing on channel.
Detector Temperature	Camera temperature out of tolerance.
Detector Locked	Camera temperature not stable.
Detector Saturation	Camera saturation < 2 % or > 80 %.
CSM Diamond Heater Temperature	Laser calibration standard temperature out of tolerance.
CSM Laser Interlock Alarm	Laser interlock interrupted, e.g. probe cable broken or improperly connected
Laser Interlock Status	Laser disabled, e.g. due to being turned off.
Laser Diode Current	Laser diode drive current approaches maximum.
Laser Cooler Open Circuit	Error received from laser.
Laser Diode Current Error	Error received from laser.
Laser Diode Temp Error	Error received from laser.
Laser Diode Temp Startup	Error received from laser.
Laser Diode Temp Warning	Error received from laser.
Laser Power Feedback	Error received from laser.
Laser Power Feedback Error	Error received from laser.
Laser Unit Temp Error	Error received from laser.
Laser Unit Temp Startup	Error received from laser.

Table 12. System warnings and errors

## 9.2.4 Quick reference troubleshooting

Diagnostic Title	Solution
Probe is not emitting laser	Check whether the probe is connected properly. Confirm that laser key is turned <b>ON</b> and indicator light is lit. Inspect whether the shutter switch is in the <b>ON</b> position. Examine whether the remote unlock connector is present on specific channel.
RunTime is frozen and not responding	Reboot device by holding the <b>ON/OFF</b> button on the front of the device for 12 seconds until it powers down. Release power button. Depress power button momentarily to power the device back up.
RunTime reports a detector temperature warning	Camera hasn't had time to cool down. The camera usually needs 20 to 25 minutes from when it powered up to cool down to the proper temperature.
The probe fiber is broken	A broken fiber cable will open the interlock and turn the laser off. For Technical Service, refer to our website ( <u>https://endress.com/contact</u> ) for the list of local sales channels in your area.
The laser has failed	Check <b>Options &gt; Diagnostics</b> for a reading on the laser's current and power. For Technical Service, refer to our website ( <u>https://endress.com/contact</u> ) for the list of local sales channels in your area.
RunTime won't initialize	Follow instructions in the Restore Console to restore an export file already saved that contains settings, calibrations, and verification data.

Table 13. Troubleshooting

## 9.3 OPC connection troubleshooting

Endress+Hauser cannot support connection to user specific OPC clients. Endress+Hauser uses a free OPC Client, UAExpert for testing server-to-client connections. A Service Engineer can provide instructions for the UA Expert Client to run Raman RunTime OPC Server tests. If problems persist, check the following items:

- Confirm whether the current version of Raman RunTime is operating.
- Verify the status of "Connected" under **Options > System > Network**.

	Netv	vork
~~~	Network 1 (CONNECTED)	00:90:FB:6F:A7:2E
503 General	DHCP	Static
ģ	IP Address	10.139.73.22 >
Analysis	Subnet Mask	255.255.255.0 >
€→	Gateway	10.139.73.1 >
utomation	Network 2 (DISCONNECTED)	00:90:F8:6F:A7:2F
吊	DHCP	Static
自由白 etwork	IP Address	>
	Subnet Mask	>
e & Time	Gateway	>
0	IDENTITY	
Security	Hostname	RAMAN-123 >
Update		
$\leftarrow$		e lu
	Ар	ріу

Figure 87. Network screen showing connected

- Make sure the correct Ethernet port is used.
- Check whether the Ethernet port is active with link/activity lights.
- Make sure a valid IP Address is shown.
- Test the ping from a direct connected laptop and ensure both are configured properly.
- Validate the OPC UA communication by using a test client such as UA Export.
- Check port configuration on client.

## 9.4 Restarting the system

In Raman RunTime, navigate to **Options > System > General** and click **Restart**. Only users with administrative privileges can restart the device.

	General	
र्ट्रे General	INSTRUMENT Instrument Name Raman	Analyzer >
Analysis	Advanced	>
$\leftarrow \rightarrow$	Calibrate Touch Screen	
Automation	Restart	
Network	Shut down	
Date & Time		
[T]		
Update		
Close	8 GB (	73% free)

Figure 88. Restart

## 9.5 Raman Rxn2 system and power loss

The instrument retains its last known power state in non-volatile memory. Should power be interrupted to the instrument at any time, the instrument will recall its last known power state and return to that power state once power is restored. For example, if the instrument was ON when power was interrupted, it will power on automatically once power is restored. If the laser was ON and the laser key is also in the ON position as well, then the laser will turn on. In the unlikely event of this occurring, this represents a possible laser exposure hazard. While power is interrupted, if the instrument was ON at the time of power loss, the power switch will flash a two-blink error code for 30 – 60 seconds, indicating that power has been lost.

## 9.6 Recovery console

The Recovery Console appears if Raman RunTime is unable to initiate. Raman RunTime will produce a prompt to enter recovery console after trying to initiate five times. There are several options for resolving the problem.



Figure 89. Recovery console menu

The Recovery Console consists of the following functions:

• Restart. This button will restart the system.



Figure 90. Restart warning

Update. This button is used to update the Raman RunTime embedded software that is provided on a USB stick.
 Put the USB flash drive in the USB slot and follow menu prompts to find the file and initiate the update.

		Select the up	date file	
	DRIVE	098-304168-USB Device (970.48 M	∕IB free space out of 14.92 G ∽	
	FOLDER	/RamanRuntimeVM-6.4.0_Beta		
	Disk			>
<				
ap on the file you wi	sh to sele	ct.		
			Cancel	
			Galifort	

Figure 91. Update mode menu

• **Service Mode.** The Service Mode option is only available to qualified service personnel. For Technical Service, refer to our website (<u>https://endress.com/contact</u>) for the list of local sales channels in your area.



Figure 92. Service mode menu

- Export. This button is used to download data, settings, or configurations of the analyzer, and only when it is in an error state. This export is intended for Endress+Hauser Support to investigate problems. In order to Restore the analyzer, a user can ONLY use a previously generated and saved (to a USB drive) Export file from within Raman RunTime. The last generated and saved export file will also reside on the system's onboard memory.
   NOTICE
- The recovery console generated export is ONLY intended for Endress+Hauser Support to investigate problems. This export file CANNOT be used to restore the analyzer.



Figure 93. Exporting a file

 Restore. This button can restore Raman RunTime to an earlier state of the analyzer that was saved at an earlier time from within Raman RunTime, and exists on the onboard storage or on a USB drive. This is the recommended solution to restoring Raman RunTime to an operational mode.

		Select the ex	port file		
	SOURCE	Onboard Exports		v	
F 2 E	laman A 022-06-2- xport	nalyzer (20220624-1903) 1 19-03-26		DATA	
arch Complete: 1	Export Fou	nd			

Figure 94. Restoring Raman RunTime to an earlier version using export stored on the analyzer storage

1. If the export file is on a USB drive, search through folder structure to find the needed file to restore settings and configuration. Please choose a Full level export file to restore the settings, configuration, models, and data. If an export is Full level and contains the aforementioned content for a full restore, this is indicated by the **Green** DATA tag next to the export file as shown below.

	Select the expor	t file	
SOURCE	098-304168-USB Device (797.9 MB free	space out of 14.92 GE 👻	
Raman /	Analyzer (20220624-1903)		DATA >
System Ex	ports		
rch Complete: 1 Export For	nd		
rch Complete: 1 Export For	nd		

Figure 95. Restoring Raman RunTime to an earlier version using export stored on a USB drive

While the settings and configuration will always be restored, Raman RunTime will prompt and ask whether the user also wants to restore data as well if a Full export (with a **Green** DATA tag) is available and selected.



Figure 96. Warning menu to select level of restore for Raman RunTime

2. Select the file to restore the analyzer. Once the file is selected, a new menu will appear to confirm the action. Click **OK**.



Figure 97. Menu to confirm restore

- 3. The system will now Restart and Restore. This can take a while, so DO NOT shut down the analyzer. Once the Restore is finished, the instrument will Restart, and Raman RunTime will Launch.
  - **Reset.** By clicking the **Reset** button, and then **Continue**, the system will reboot and delete all of the calibrations, models, acquisitions, logs, and other data.



*Figure 98. Menu to confirm resetting the device*
# 10 Maintenance

## 10.1 Optimization

If you move the Raman Rxn2, it may be necessary to re-optimize its performance. First, re-verify its performance using Raman RunTime and compare the results for the current to the previous verification. If the signal intensity has dropped significantly, you may benefit from following the optimization guidelines below.

### 10.1.1 Sample position

If the sample has been moved from the focal point of the probehead, less Raman scattering is recovered by the probehead lens and transmitted to the spectrograph. This is the easiest area to check first.

Perform the following procedure in a darkened room:

- 1. Click Focus on the Stream Detail View.
- 2. Watch as the signal increases and decreases in response to the sample movement in front of the probehead lens.
- 3. Be aware of any potential laser light reflected off the sample container during this procedure.

### 10.1.2 Cleaning the lens or window

If the lens or window on the probehead has come in contact with a sample, dust or fingerprints, etc., it may need to be cleaned. Use a lens cloth and a water-based lens-cleaning solution to gently wipe away any contaminant.

### 10.1.3 Detector camera alignment

If the Raman Rxn2 spectrograph internal optics have shifted, the detector camera alignment may need to be changed.

#### 

• The CCD camera alignment is set at the factory and rarely needs to be altered in the field. Alignment should be performed only by experienced personnel.

Prior to performing a camera alignment operation, it is important to ensure that no stray light is entering any of the probes that are attached to the Raman Rxn2. Alignment is performed with an internal white light source, and stray light entering on any of the attached probes can interfere with the alignment light source.

To perform the camera alignment:

- 1. Navigate to **Options > Calibration**.
- 2. Click **Calibrate** under the Internal Calibration section then select **Recalibrate All** from the Calibration Mode drop-down list. Click **Calibrate**.

#### NOTICE

• All probe calibrations and verifications are invalidated after Recalibrate All and will need to be re-performed.

#### 10.1.4 Calibration and verification wizard

Reliable, transferable calibration is important for comparing data acquired at various times or with different analyzers. Different instruments analyzing the same sample can generate nearly identical spectra if they are properly calibrated. The Raman RunTime software package includes an automatic calibration wizard that guides you through a procedure to automatically calibrate the wavelength and intensity axes and the wavelength of the laser.

#### NOTICE

• After the initial calibration during installation, the Calibrate Periodically function is usually sufficient to maintain the wavelength and laser calibration of the Raman Rxn2.

A summary of the recommended calibration and verification sequence is as follows:

- 1. Internal Calibration.
- 2. Probe Calibration (Raman Calibration Accessory).
- 3. Probe Verification (external 70 % IPA or cyclohexane sample where applicable).

#### 10.1.4.1 Internal calibration

Internal calibration includes spectrograph alignment, spectrograph x-axis wavelength, and laser wavelength calibrations.

The wavelength calibration of the Raman Rxn2 can be automatically checked at a frequency of choice by navigating to **Options > Calibration > Internal Calibration** and selecting **Calibrate Periodically**. At least once per day is recommended. Once enabled, the Raman Rxn2 will automatically compare the current analyzer response (spectrograph and laser wavelength calibrations) to calibration specifications and will recalibrate if the spectrograph wavelength or laser wavelength is out of spec.

#### 10.1.4.2 Probe calibration

The sensitivity of the Raman Rxn2 varies with wavelength due to variations in the throughput of the optics and the quantum efficiency of the CCD. The Probe Calibration function in your embedded software application can be used to remove the effects of this variation from measured spectra. Probe calibration can be performed using the automatic calibration wizard in Probe Calibration on the Raman Rxn2 requires the Raman Calibration Accessory.

Attach the lamp head to the probe on channel 1 using the appropriate optic adapter. Each channel must be probecalibrated separately. Follow the instructions in the automatic calibration wizard.

#### 10.1.4.3 Probe verification

#### 

Bio probes such as bIO-Optic, Rxn-45, and single-use adapters, should ONLY use 70 % IPA. Only 70 percent by volume (%v/v) will work. Endress+Hauser recommends using CiDehol 70 by Decon Laboratories. Other probes can use cyclohexane.

The Probe Verification wizard may be used to verify that the Raman Rxn2 is performing within specifications. Probe verification will acquire a Raman spectrum of a standard Raman sample, usually 70 % IPA or cyclohexane, and analyze the resultant spectrum for peak positions, peak area ratios and Raman signal strength. The peak position verification will verify that the spectrograph and laser wavelength calibrations are within specification. The peak area ratio verification will verify that the probe intensity calibrations are within specification. The signal strength verification will verify that the instrument's signal-to-noise ratio is within specification. A report will be generated showing the results of the verification steps along with a Pass/Fail indication.

## 10.2 Servicing the Raman Rxn2 analyzer

This section is divided into two parts: 1) embedded software-related symptoms and 2) hardware-related symptoms. The distinction is not clear-cut in many cases, so check the other section if your situation is not described in the expected section.

Some of the service procedures require removal of the protective cover. Special precautions are therefore necessary to deal with the added optical and electrical hazards that are present during service operations.

#### 

General users should not open the Raman Rxn2 enclosure. Only qualified personnel familiar with high voltage electronics should open the system enclosure to perform necessary maintenance or service. Opening the Raman Rxn2 enclosure can expose users to hazardous laser radiation or high voltage electronics.

#### 10.2.1 List of embedded software related symptoms

#### Symptom: Raman spectra contain non-random noise

**Possible Solutions:** 

1. The intensity calibration file is no longer valid.

The intensity calibration file is a map of the total instrument response of the system (CCD quantum efficiency, grating and lens efficiencies, etc.). If the spectrograph focal plane shifts with respect to the CCD detector, then the intensity calibration map is no longer correct. An incorrect intensity calibration file will not remove pixel-to-pixel sensitivity variation, which is as high as 3 % for some CCD chips. Shifting of the spectrograph focal plane on the CCD detector can result from adjusting the Raman Rxn2 spectrograph alignment, mechanical shock applied to the Raman Rxn2 or substantial ambient temperature shifts. Recalibrate the intensity axis.

2. The intensity per pixel per accumulation exceeded the half-well potential of CCD pixels during the creation of the intensity calibration file.

Non-linearity reduces the effectiveness of the intensity calibration file at correcting the CCD pixel-to-pixel sensitivity variation. Recalibrate the intensity axis.

#### Symptom: Intensities in Raman spectra show non-linear behavior or Raman peak shapes are distorted

**Possible Solutions:** 

1. Repeat the data acquisition using a shorter data acquisition time and perhaps more accumulations (which adds intensities in the embedded software rather than on the CCD chip).

The signal from the CCD can become slightly non-linear as the photo-generated charge approaches the full-well potential. This can cause non-linear behavior and peak distortion in Raman spectra.

2. The intensity calibration file is invalid.

If the intensity calibration was made using a light source that did not evenly illuminate the entire collection lens of the spectrograph, then any experiment that does not reproduce the exact uneven illumination at the time of the intensity calibration will not have the same instrument response and therefore will not be properly corrected. Recalibrate the intensity axis.

#### Symptom: Raman RunTime reports a detector temperature warning

**Possible Solutions:** 

1. The camera has not had time to cool down.

The camera usually requires 15 to 20 minutes from the time it is powered up until it has cooled down to the proper temperature.

### 10.2.2 List of hardware-related symptoms

#### Symptom: Several sharp artifact peaks appear in all Raman spectra

Possible Solutions:

1. The room lights are introducing an emission line spectrum into the spectrograph.

Turn off any fluorescent lights during future experiments. Cover reactor with aluminum or other light-blocking material to prevent light entering.

#### Symptom: The number of intensity counts from verification is significantly less than specification

Possible Solutions:

1. The sample is not in the focal plane of the fiber optic probehead.

Adjust the position of the sample with respect to the probe focus.

- 2. The fiber optic cable is not attached correctly to the Raman Rxn2 base unit. Check fibers for proper engagement and that they are latched.
- 3. The laser power reaching the sample is too low.

Measure the laser power at the sample and compare it to the normal power for your configuration. For Technical Service, refer to our website (<u>https://endress.com/contact</u>) for the list of local sales channels in your area.

#### Symptom: Spectrum sits on top of a broad featureless halo

Possible Solutions:

1. Probe tip may be fouled.

Pull probe from process and clean the probe per the appropriate service manual instructions. For Technical Service, refer to our website (<u>https://endress.com/contact</u>) for the list of local sales channels in your area.

#### Symptom: No laser light reaches the sample

**Possible Solutions:** 

1. The laser is not on.

Check to see if the laser key is in place and the indicator is lit.

2. The laser is not lasing.

For Technical Service, refer to our website (<u>https://endress.com/contact</u>) for the list of local sales channels in your area.

3. The internal power plug has disconnected from the laser.

Open the Raman Rxn2 base unit enclosure. The computer-style power cord should be firmly plugged into its socket on the laser.

4. The fiber optic cable is not attached correctly to the Raman Rxn2 base unit.

Check hybrid fibers for proper engagement and that they are latched.

5. The laser injector is misaligned.

For Technical Service, refer to our website (<u>https://endress.com/contact</u>) for the list of local sales channels in your area.

6. The remote interlock connector shorting plug is unplugged.

Make sure the remote interlock connector shorting plugs for all channels are installed. Make sure the corresponding front panel interlock indicators are illuminated.

7. The probe fiber is broken.

The interlock cuts power to the system if the cable is broken. If this happens, contact Endress+Hauser or an authorized distributor.

8. The Calibration Switching Module has failed.

For Technical Service, refer to our website (<u>https://endress.com/contact</u>) for the list of local sales channels in your area.

#### 9. The laser has failed.

Check **Options > Diagnostics** for laser diode current and power.

#### Symptom: Embedded software indicates an excessive number of tracks found during recalibrate all

**Possible Solutions:** 

1. Stray light from connected probes are entering spectrograph during camera alignment.

Cover all connected probe optics to prevent any stray light from entering the spectrograph.

#### Symptom: Embedded software indicates a failure during wavelength calibration

**Possible Solutions:** 

1. Stray light from connected probes is entering spectrograph during calibration.

Cover all unused but connected probe optics to prevent any stray light from entering the spectrograph. Also, ensure that probes being used for sampling are protected from stray light. Wavelength calibration is performed with a light source internal to the Raman Rxn2 base unit. If stray light from connected probes can enter the spectrograph, it may interfere with the internal calibration lamp.

#### Symptom: Embedded software indicates internal temperature error

- 1. The filter needs attention.
- Clean or replace the filter.
- 2. The ambient temperature is above 35°C (95 °F).

Lower ambient temperature to a temperature within specified ambient temperature range.

## 10.3 Support

Diagnostic exports and the current Raman RunTime version number will be needed for Endress+Hauser Support. The version can be viewed from **Options > System > Update**.

Current Version: 6.4.(11B2D0E0+)

Click Information to display additional details about the versions and updates applied.

	Current Version:	6.4 (41B2D9E9+)	
Runtime Core Profile:	e: 6.4 (41B2D9E9+) 785 Rxn		Â
Jpdates Inst	alled:		
ResetDevice			
2022-06-20 13	:37	ОК	
JpdateSSLCe	rtificates		
2022-06-20 07	:14	ОК	
JpdateOpcUa	Certificates		
2022-06-20 07	:13	ОК	
SyncTime			
2022-06-20 07	:13	ОК	
JpdateSSLCe	rtificates		
2022-06-20 07	:08	ОК	
JpdateOpcUa	Certificates		
2022-06-20 07	:07	ОК	
SyncTime			_

Figure 99. Current version

### **10.3.1** Contact information

For Technical Service, refer to our website (<u>https://endress.com/contact</u>) for the list of local sales channels in your area.

# 11 Repair

## 11.1 Servicing and spare parts

The two main user-serviceable or replaceable parts are the disposable air filter and the laser. Part numbers for these items are listed in the table below. The laser assembly has been designed for easy replacement and this is typically accomplished by the customer. Optionally, an Endress+Hauser service engineer can install the laser during a contracted service visit.

#### NOTICE

• The performance of procedures (including service), the use of controls, or the adjusting of the instrument other than as specified in the manual will void the warranty.

Model Number	Description	
RXN2 Air Filter	One packet of replacement Air Filters for a Raman Rxn2 analyzer (Quantity 5 filters)	
ULA-EM-785-2/4	Embedded NIR 785 nm Diode Laser for Raman Rxn2 features:	
	<ul> <li>Laser Wavelength - 785 nm</li> <li>&gt;125 mW of 785 nm laser power delivered to the Probehead*</li> <li>Integrated Holographic Laser Bandpass Filter</li> <li>Universal Laser Injector Assembly</li> <li>One Year Unlimited Hours Warranty</li> </ul>	
	*Using standard multi-mode optical fiber	
L-INV-EM2/4-785-R	Refurbished Embedded NIR 785 nm Diode Laser for Raman Rxn2 analyzer features:	
	<ul> <li>Return of Existing NIR 785 nm Diode Laser and Injector Assembly</li> <li>Clean and Refurbish Laser and Injector as necessary</li> <li>Refurbished Specifications:         <ul> <li>&gt;150 mW of 785 nm laser power delivered to the Probehead*</li> <li>One Year Unlimited Hours Laser Warranty</li> </ul> </li> </ul>	
	*Using standard multi-mode optical fiber	
	Requires Trade-in of Existing Customer-owned Embedded NIR 785 nm Diode Laser (serial number to be provided)	
ULA-EM-532-2/4	Embedded 532 nm Frequency Doubled Diode-Pumped Nd:YAG Laser for Raman Rxn2 analyzer features:	
	<ul> <li>Laser Wavelength - 532 nm</li> <li>Diode-Pumped Nd:YAG Laser</li> <li>150 mW Laser Head Output</li> <li>One Year / 5000 Hours Warranty</li> <li>&gt;80 mW of 532 nm laser power delivered to the probehead*</li> </ul>	
	*Using standard multi-mode optical fiber	
L-INV-EM2/4-532/K	<ul> <li>Refurbished Embedded VIS 532 nm Diode Laser for Raman Rxn2 features:</li> <li>Return of Existing VIS 532 nm Diode-Pumped Nd:YAG Laser and Injector Assembly</li> <li>Clean and Refurbish Laser and Injector as necessary</li> <li>Refurbished Specifications: <ul> <li>&gt;80 mW of 532 nm laser power delivered to the probehead*</li> <li>One Year or 5000 Hours Laser Warranty</li> </ul> </li> </ul>	
	*Using standard multi-mode optical fiber	
	Requires Trade-in of Existing Customer-owned Embedded VIS 532 nm Diode-Pumped Nd:YAG Laser (serial number to be provided)	
ULA-EM-1000-2/4	Embedded NIR 993 nm Diode Laser for Raman Rxn2 analyzer features:	
	<ul> <li>Laser Wavelength – 1000 nm</li> </ul>	

The following table provides a list of common parts that may be ordered and installed.

N# 1 1 NY 1		
Model Number	Description	
	<ul> <li>&gt;150 mW of 1000 nm laser power delivered to the Probehead*</li> </ul>	
	<ul> <li>Integrated Holographic Laser Bandpass Filter</li> </ul>	
	<ul> <li>Universal Laser Injector Assembly</li> </ul>	
	<ul> <li>One Year Unlimited Hours Warranty</li> </ul>	
	*Using standard multi-mode optical fiber	
L-INV-EM2/4-1000/R	Refurbished Embedded NIR 1000 nm Diode Laser for Raman Rxn2 analyzer features:	
	<ul> <li>Return of Existing NIR 1000 nm Diode Laser and Injector Assembly</li> </ul>	
	<ul> <li>Clean and Refurbish Laser and Injector as necessary</li> </ul>	
	<ul> <li>Refurbished Specifications:</li> </ul>	
	<ul> <li>&gt;150 mW of 1000 nm laser power delivered to the Probehead*</li> </ul>	
	One Year Unlimited Hours Laser Warranty	
	*Using standard multi-mode optical fiber	
	Requires Trade-in of Existing Customer-owned Embedded 1000 nm Laser (serial number to be	
	provided)	

Table 14. Spare parts

For Technical Service, refer to our website (<u>https://endress.com/contact</u>) for the list of local sales channels in your area.

# 12 Technical data

## 12.1 Specifications

Raman Rxn2 analyzers may be configured to operate with one of several different laser wavelengths. Currently, Raman Rxn2 analyzers may be equipped with a 532 nm, 785 nm, or 993 nm laser as standard.

## 12.1.1 Base unit

Item	Description
Operating temperature	15 °C to 30 °C (59 °F to 86 °F)
Storage temperature	–15 °C to 50 °C (5 °F to 122 °F)
Relative humidity	20–80 %, non-condensing
Warm-up time	120 minutes
Operating voltage	100-240 V, 50-60 Hz, ±10 %
Transient over-voltages	Over-voltage category 2
Power consumption (maximum)	400 W
(typical startup)	250 W
(typical running)	120 W
Base Unit Dimensions (width x depth x height)	279 × 592 × 483 mm (11 in. × 23.3 in. × 19 in.)
Optional Cart Dimensions	71 cm × 76 cm x 101.5 cm to tabletop
(width x depth x height)	147 cm to monitor top
	(28 in. wide × 30 in. deep x 40 in. to tabletop)
	(58 in. to monitor top)
Weight	
Base Unit	71 lbs. (32 kg)
Cart Mounted	205 lbs. (93 kg)
Sound Level (from operator's perspective)	58.9 dB, A-weighted
IEC 60529 rating	IP20
Altitude	Up to 2000m
Pollution degree	2

Table 15. Base unit

## 12.1.2 Spectrograph

Item	Description
Туре	Proprietary axial transmissive
Aperture ratio	<i>f</i> /1.8
Focal length	85 mm
Grating (1 or 4 channels, 532 nm, 785 nm)	HoloPlex transmissive (Rxn2-785 hybrid analyzers use HoloSpec transmissive)
Grating (1 or 4 channels, 1000 nm)	HoloSpec transmissive
Spectral coverage (532 nm)	150 to 4350 cm <sup>-1</sup>
Spectral coverage (785 nm)	150 to 3425 cm <sup>-1</sup>
Spectral coverage – Rxn2 hybrid (785 nm)	175 to 1890 cm <sup>-1</sup>
Spectral coverage (1000 nm)	200 to 2400 cm <sup>-1</sup>
Slit	50 µm fixed
Spectral resolution (532 nm)	5 cm <sup>-1</sup> average
Spectral resolution (785 nm)	4 cm <sup>-1</sup> average
Spectral resolution (1000 nm)	5 cm <sup>-1</sup> average

Table 16. Spectrograph

## 12.1.3 Laser

Item	Description
532 nm	
Excitation wavelength	532 nm
Maximum power output	120 mW
Warranty	unlimited hours for 1 year
785 nm	
Excitation wavelength	785 nm
Maximum power output	400 mW
Warranty	unlimited hours for 1 year
1000 nm	
Excitation wavelength	993 nm
Maximum power output	400 mW
Warranty	unlimited hours for 1 year

Table 17. Laser

## 12.1.4 Probes

Item	Description
Raman Rxn2 (single channel, four channel,	Compatible with:
and starter)	Rxn-10 probehead equipped with immersion or non-contact optics
	Endress+Hauser gas phase Raman probe
	Endress+Hauser liquid phase Raman probes
	Endress+Hauser bioprocessing Raman probes
Raman Rxn2 hybrid	Compatible with:
	Rxn-20 probe and
	One other ALT probe including:
	Rxn-10 probehead equipped with immersion or non-contact optics
	Endress+Hauser gas phase Raman probe
	Endress+Hauser liquid phase Raman probes
	Endress+Hauser bioprocessing Raman probes

## 12.2 Hazard area certification

Endress+Hauser analyzers carry IECEx, ATEX, and NA certifications for installation in general purpose zone with output into hazardous areas. For more specific information on the hazardous area rating for field measurements, refer to the operating manual of the installed probe.

Item	Description
Certification – Base Unit (fiber optic and interlock outputs only)	ATEX         Image: Control (Control (Contro) (Contro) (Control (Control (Control (Control (Contro
	IECEx ITS 20.0050X
	<b>North American</b> Intrinsically Safe output for probe and sensors used in Class I, Division 1, Groups A, B, C, and D. Ambient Temperature Range: $5^{\circ}$ C (41 °F) $\leq$ Tamb $\leq$ 35°C (95 °F)

Table 18. Hazardous area certification

# **13** Supplementary documentation

All documentation is available:

- On the media device supplied (not included in the delivery for all device versions).
- Smart phone/tablet: Endress+Hauser Operations App.
- For Technical Documentation, refer to our website (<u>https://endress.com/downloads</u>).

Part Number	Document Type	Document Title
KA01544C	Brief Operating Instructions	Raman Rxn2 Brief Operating Instructions
XA02700C	Safety Instructions	Raman Rxn2 Safety Instructions
TI01608C	Technical Information	Raman Rxn2 Technical Information

Table 19. Supplementary documentation

# 14 Index

ac power, 32 safety, 10 air exhaust, 28 filter, 39, 78 flow, 19 temprature, 19 ventilation, 19 altitude, 80 analyzer calibration window, 26 four channel, 12 hybrid, 12 one channel, 12 probe, 20 rear panel, 16 status, 66 aperture laser, 38 numeric, 10 ratio. 81 asynchronous updates, 30 automatic calibration, 41 calibration wizard, 73 connection, 29 dark exposures, 47 user privilege, 60 automation, 29 base unit air filter, 39 enclosure, 19 humidity, 19 laser, 38 lifting, 8 raman rxn2, 12 rear panel, 16 specifications, 80 temperature, 19 ventilation, 19 basic connectivity, 29 batch data export, 56 calibration internal, 41, 74 module, 38 probe, 74 probes, 42 reports, 46 verification, 26 wizard, 73 camera alignment, 73

interface, 67 saturation. 68 temperature, 68 cdrh compliance, 11 certification compliance label, 11 hazardous area, 82 ua. 58 clean, 73 commissioning, 29 compliance labels, 11 conditions of use, 11 continuous mode, 51 dark exposures, 47 data batch export, 56 collection, 46 tag, 72 technical, 80 transfer, 56 dc fuses. 38 safety, 10 diagnostics environment, 63 exports, 65 trends, 64 dimensions, 80 electrical block diagram, 33 connection, 28 safety, 10 errors, 66 export, 62 batch data, 56 compliance, 4 diagnostics, 65 spc file network, 56 external, 31 eye optical safety, 10 protection, 9 focal length, 81 focus. 48. 73 four channel analyzer, 12 calibration window, 26 probe, 20 rear panel, 16 front panel, 14 fuses, 38 glossary, 5 grating, 81

hca. 19 health. 9 humidity, 19 hybrid analyzer, 12 calibration window, 26 probe, 20 rear panel, 16 i/o panel, 31 interlock connector, 11, 36 key control, 11 laser apetures, 38 base unit, 38 interlock circuit, 38 key switch. 14 low power, 67 safety, 9 safety notice, 9 lens, 73 lifting raman rxn2, 8 manual mode, 50 mini display port, 31 mobile cart, 39 modbus, 58 models loading, 53 results. 54 network configuration, 25 numeric aperture, 10 on/off button flashing, 67 opc connection, 69 ua. 29 operating voltage, 80 optical safety, 10 optimization, 73 password management, 61 periodic mode, 52 pollution degree, 80 port ethernet, 69 exc. 24 rs-485, 16 usb, 14 power consumption, 80 electrical, 19 grounding, 32 switch, 14 up, 25 probe calibration, 42 connection status, 14 laser apetures, 38 on/off, 9 verification, 74

protective housing, 11 raman runtime initial setup, 40 network configuration, 25 overview, 13 raman rxn2 air filter, 39 humidity, 19 interior, 37 lifting, 8 location, 19 temperature, 19 ventilation, 19 rear panel, 16 recovery console, 70 relative humidity, 19 remote interlock connector, 11, 36 remote access, 56 repair, 78 restart, 69 safety electrical. 10 health, 9 optical, 10 sample position, 73 security additional features, 61 basic connectivity, 29 password, 61 software, 59 user privileges, 60 single channel analyzer, 12 calibration window, 26 probe, 20 rear panel, 16 slit, 81 snapshot mode, 49 software raman runtime, 13, 40 requirements, 62 security, 59 user privilege, 60 sound level, 80 spare parts, 78 spc file network, 56 hvbrid. 44 probe calibration, 42 specification, 80 spectrograph, 38 status bar, 66 status indicator, 57 support, 77 symbols, 4 temperature, 19

test server, 69 touchscreen, 31 trends, 64 us export compliance, 4 usb port, 14 user privileges, 60 ventilation, 19 verification calibration, 26 probe, 74 reports, 46 wizard, 73 warm-up time, 80 warnings, 66 weee directive, 11 weight, 80

www.addresses.endress.com



People for Process Automation