



















# FMP 40 in Level Measurement in Bioreactor-Pharm.

# Levelflex M FMP 40 in pharmaceutical reactor vessels



Typical small SS tanks

The guided wave radar Levelflex M FMP 40 is used to measure level in a bioreactor.

## Customer profile

A large producer of pharmaceuticals and biologics in California.

## Application description

Product: Various enzymes, WFI Temperature: Ambient

Stainless steel, 40" tall Tanks:

Tri-clamp Mounting: Dielectric constant: >10



FMP40 with TSP Tri-clamp fitting

# Application challenge

The customer used to use magneto strictive transmitters but had problems with the accuracy due to changing densities of the product and its cleanability. High accuracy is required, plus the customer wanted to calibrate the instrument and be able to change electronics without opening the tank.

# Application solution

A Levelflex M FMP 40 with Tri-clamp was installed in the tank. The measurement results showed excellent measuring, even with changing densities and foam in the tank. The special bio-design of the transmitter made it easy to remove the electronics from the probe, place it on a reference probe and calibrate without breaking the sterility of the reactor. Fluctuations in density, temperature and foam do not affect the measurement.



Checking levels in bio-storage vats

#### Instrument description

The Levelflex M is a top mounted compact level transmitter that transmits micro impulse radar signals on a rod or cable. The pulses are reflected by the product surface, received by the electronics, evaluated and converted into level information.

Wetted materials are FDA, USP 88 Class VI. The rod is electropolished and the seal

For more information, contact Endress+Hauser, Inc. 317-535-7138 www.us.endress.com





FMP 40 with Tri-clamp allows electronics to be removed without disturbing process.



FMP 40 Tri-clamp assembly

ISO 9001:2000 Certified

# USA

Endress+Hauser, Inc. 2350 Endress Place Greenwood, IN 46143 Tel. 317-535-7138 Sales 888-ENDRESS Service 800-642-8737 Fax 317-535-8498 inquiry@us.endress.com www.us.endress.com

#### Canada

Endress+Hauser, Canada 1075 Sutton Drive Burlington, ON L7L 5Z8 Tel. 905-681-9292 800-668-3199 Fax 905-681-9444 info@ca.endress.com www.ca.endress.com

## Mexico

Endress+Hauser México, S.A. de C.V.
Fernando Montes de Oca 21 Edificio A Piso 3
Fracc. Industrial San Nicolas
54030. TlaInepantla de Baz
Estado de México
México
Tel. +52 55 5321 2080
Fax +52 55 5321 2099
eh.mexico@mx.endress.com
www.mx.endress

